

Finnigan[™]
ConFlo III
Universal Interface

Operating Manual

Revision B
111 0042

*Finnigan, Flash EATM and Eager 300 are trademarks of Thermo Electron Corporation.
CE Instruments and Carlo Erba Instruments are part of Thermo Electron Corporation.
Thermo Electron (Bremen) and Thermo Electron (Italy) are part of Thermo Electron Corporation.
Swagelok[®] is a registered trademark of the Crawford Fitting Company.
Teflon[®] is a registered trademark of E. I. du Pont de Nemours & Co.
Windows 95, Windows 98, Windows 2000, Windows NT and Windows XP are trademarks of Microsoft Corporation.*

Technical information contained in this publication is for reference purposes only and is subject to change without notice. Every effort has been made to supply complete and accurate information; however, Thermo Electron assumes no responsibility and will not be liable for any errors, omissions, damage, or loss that might result from any use of this manual or the information contained therein (even if this information is properly followed and problems still arise).

This publication is not part of the Agreement of Sale between Thermo Electron and the purchaser of a Thermo Electron system. In the event of any conflict between the provisions of this document and those contained in Thermo Electron Terms and Conditions, the provisions of the Terms and Conditions shall govern.

Reference to System Configurations and Specifications supersede all previous information and are subject to change without notice.

Printing History: Revision A printed as Draft Version in August 2003.

Revision B printed in December 2003.

Thermo

ELECTRON CORPORATION

The products of Thermo Electron Bremen are produced under ISO 9001 accredited quality management systems.

Australia: P.O. Box 239 Rydalmere • Unit 14, 38 – 46 South Street • Rydalmere, N.S.W. 2116 • [61] (02) 9898-9000
Austria: Wehlstrasse 27b • A-1200 Wien • [43] (01) 333 50 34-0
Belgium: Technologiestraat 47 • B-1082 Brussels • [32] (02) 482 30 30
Canada: 5716 Coopers Avenue, Unit 1 • Mississauga, Ontario • L4Z 2E8 • [1] (905) 712-2258
France: 16 Avenue du Québec • Silic 765 • Z.A. de Courtaboeuf • F-91963 Les Ulis Cédex • [33] (01) 60 92 48 00
Germany: Im Steingrund 4-6 • D-63303 Dreieich • [49] (06103) 408 0
Italy: Strada Rivoltana • I-20090 Rodano (Milano) • [39] (02) 95059 226
Japan: C-2F • 3-9, Moriya-cho, Kanagawa-ku • Yokohama, Kanagawa • 221-0022 • [81] (45) 453 9100
Japan: Esaka Grand Building • 2-3-1 Esaka-cho, Suita City • Osaka 564-0063 • [81] (06) 6387-6681
Netherlands: Takkebijsters 1 • 4817 BL Breda • [31] (076) 5878 722
P.R. China: Room 901, Ping-an Mansion • No. 23, Jin Rong Street • Xi Cheng District • Beijing 100032 • [86] (010) 6621 0839
Spain: Sepulveda 7 A • ES-28108 Alcobendas (Madrid) • [34] (091) 657 4930
Spain: Acer 30 – 32 • Edificio Sertram – Planta 2, Modulo 3 • ES-08038 Barcelona • [34] (093) 223 0918
Sweden: Pyramidbacken 3 • S-141 75 Kungens Kurva (Stockholm) • [46] (08) 556 468 00
United Kingdom: Stafford House • 1 Boundary Park • Boundary Way • Hemel Hempstead • Hertfordshire HP2 7GE • [44] (01442) 233 555
U.S.A.: 355 River Oaks Parkway • San Jose, CA 95134-1991 • [1] (408) 965-6000

Notes: The country code is enclosed in square brackets []. The city code or area code is enclosed in parenthesis (). For countries other than the U.S.A., when you are dialing from within the specified country, dial the 0 of the city code. For countries other than Italy, when you are dialing from outside the country, do not dial the 0 of the city code.

Thermo

ELECTRON CORPORATION

Customer Registration... Register now and receive all the privileges associated with being a Thermo Electron, Finnigan product user, including application reports and technical reports.

Name _____

Title _____

Company _____

Address _____

City/State/Postal Code _____

Country _____

Telephone _____ **Ext.** _____

Tell us more... Let us know more about how you use this product:

My Organization Is: (Check one only)

- Commercial (for profit) lab
- Government lab
- Hospital / Clinic
- Research Institute
- University / College
- Veterinary
- Other _____

Job Function: (Check one only)

- Administration
- Lab Management
- Operator
- Other _____

My Primary Application Is: (Check one only)

- Analytical
- Biomedical
- Clinical / Toxicology
- Energy
- Food / Agriculture
- Forensic / Toxicology
- Pharmaceutical
- Research / Education
- Other _____

Reader Survey... Help us to improve the quality of our documentation by answering a few questions:

*ConFlo III Universal Interface
Operating Manual*

*Revision B
111 0042*

	Strongly Agree	Agree	Disagree	Strongly Disagree
The manual is well organized.	1	2	3	4
The manual is clearly written.	1	2	3	4
The manual contains all of the information I need.	1	2	3	4
The instructions are easy to follow.	1	2	3	4
The instructions are complete.	1	2	3	4
The technical information is easy to understand.	1	2	3	4
The figures are helpful.	1	2	3	4

Additional Comments: (Attach additional sheets if necessary.)

Tear this sheet from the manual, fold it closed, stamp it, and drop it in the mail.

From _____

Place
Stamp
Here

**Thermo Electron (Bremen) GmbH
Finnigan Advanced Mass Spectrometry
Product Marketing
Barkhausenstr. 2**

**D-28197 Bremen
Germany**

← fold

← fold

Reparatur - Begleitkarte

Bitte vollständig ausfüllen

Repair-Covering Letter

Please fill in completely

Absender:
Dispatcher:

Geräte-Type:
Instrument Type:

Service-Nr.:
Service-No.:

Sie erhalten zur Reparatur unter unserer Bestell-Nr.:
You receive for repair under our Order No.: _____

- DOA**
- Teil wurde nicht benutzt / part not used
- Teil wurde nur für Testzwecke verwendet / used for test purposes only
- Festgestellte Mängel oder deren Auswirkung:
Established Defect or its Effect:**

Bitte detaillierte Angaben machen / Please specify in Detail

Ein Austauschteil haben wir erhalten unter Kommissions-Nr.:
An Exchange Part already received with Commission No.:

Ja/Yes Nein/No

Die Anlage ist außer Funktion
The System is out of Function

Ja/Yes Nein/No

Durch die nachfolgende Unterschrift bestätige(n) ich/wir, daß die o.g. Teile frei von gesundheits-
schädlichen Stoffen sind, bzw. vor Ihrer Einsen-
dung an Thermo Electron Bremen dekontami-
niert wurden, falls die Teile mit giftigen Stoffen in
Verbindung gekommen sind.

By signing this Document I am/we are certifying
that the a.m. Parts are free from hazardous
materials. In case the parts have been used for
the analysis of hazardous substances I/we
attest that the parts have been decontaminated
before sending them to Thermo Electron
Bremen.

Datum / Date

Unterschrift / Signature

**Notice on the Proper Use of
Thermo Electron Bremen Instruments**

In compliance with international regulations: If this instrument is used in a manner not specified by Thermo Electron Bremen, the protection provided by the instrument could be impaired.

Safety and EMC Information

In accordance with our commitment to customer service and safety, these instruments have satisfied the requirements for the European CE Mark including the Low Voltage Directive.

Designed, manufactured and tested in an ISO9001 registered facility, this instrument has been shipped to you from our manufacturing facility in a safe condition.

Caution. This instrument must be used as described in this manual. Any use of this instrument in a manner other than described here may result in instrument damage and/or operator injury.

Identifying Safety Information

This reference manual contains precautionary statements that can prevent personal injury, instrument damage, and loss of data if properly followed. All statements of this nature are called to your attention through the use of bold type and the following icons:



Warning!



Cold Surface!



Hot Surface



Strong
Magnetic Field!



High Voltage!

Every instrument has specific hazards, so be sure to read and comply with the following precautions. They will help ensure the safe, long-term use of your system.

1. Before plugging in any of the instrument modules or turning on the power, always make sure that the voltage and fuses are set appropriately for your local line voltage.
2. Only use fuses of the type and current rating specified. Do not use repaired fuses and do not short-circuit the fuse holder.
3. The supplied power cord must be inserted into a power outlet with a protective earth contact (ground). When using an extension cord, make sure that the cord also has an earth contact.

4. Do not change the external or internal grounding connections. Tampering with or disconnecting these connections could endanger you and/or damage the system.

Caution. The instrument is properly grounded in accordance with regulations when shipped. You do not need to make any changes to the electrical connections or to the instrument's chassis to ensure safe operation.

5. Never run the system without the housing on. Permanent damage can occur.
6. Do not turn the instrument on if you suspect that it has incurred any kind of electrical damage. Instead, disconnect the power cord and contact a Service Representative for a product evaluation. Do not attempt to use the instrument until it has been evaluated. (Electrical damage may have occurred if the system shows visible signs of damage, or has been transported under severe stress.)
7. Damage can also result if the instrument is stored for prolonged periods under unfavorable conditions (e.g., subjected to heat, water, etc.).
8. Always disconnect the power cord before attempting any type of maintenance.
9. Capacitors inside the instrument may still be charged even if the instrument is turned off.
10. The superconducting magnet is still charged even if the instrument is turned off.
11. Never try to repair or replace any component of the system that is not described in this manual without the assistance of your service representative.

Contents

Read This First	v
Changes to the Manual and Online Help	vi
Abbreviations	vii
Typographical Conventions	xi
Data Input	xi
Boxed Information.....	xii
Topic Headings.....	xiii
Reply Cards	xiv
Prior to Measurement	1-1
1.1 Precautions.....	1-2
1.2 About this Manual.....	1-3
1.3 Introduction.....	1-4
The ConFlo III Interface	2-1
2.1 Introduction - The Open Split Principle.....	2-2
Procedure.....	2-2
The Sample Section.....	2-3
The Reference Section.....	2-4
2.2 Front and Rear Panel of ConFlo III	2-6
Front Panel.....	2-6
Side Views	2-7
Rear Panel and Gas Connections.....	2-8
Gas Flow Diagram.....	2-9
Spare Parts of Reference and Sample Sections	2-10
Manometer.....	2-11
ConFlo III Interface and Elemental Analyzer	3-1
3.1 ConFlo III Software in Old Isodat	3-2
3.2 ConFlo III Configuration in Isodat 2.0	3-3
3.3 Installation of the A200S and GC PAL Autosampler	3-4
3.4 Manual Operation of ConFlo III.....	3-8

3.5	How to Get Started with ConFlo III and Elemental Analyzer	3-10
Nitrogen Measurement		4-1
4.1	Zero Enrichment of Nitrogen (Standard On/Off Test).....	4-2
	Instrument	4-2
	Time Events	4-3
	Evaluation	4-4
	Peak Detection	4-5
	Printout.....	4-6
4.2	Linearity Test of Nitrogen	4-9
Carbon Measurement		5-1
5.1	Zero Enrichment of Carbon (Standard On/Off Test).....	5-2
	Instrument	5-2
	Time Events	5-3
	Evaluation	5-4
	Peak Detection	5-5
	Printout.....	5-6
5.2	Linearity Test of Carbon.....	5-9
Dual Measurement		6-1
6.1	Introduction	6-2
6.2	Procedure.....	6-3
6.3	How to Perform a Jump Calibration.....	6-4
	Jump Calibration - Before Start	6-6
	Comment: If no Jump Calibration is Available	6-7
	Jump Calibration - Procedure	6-9
Sulfur Measurement		7-1
7.1	Introduction	7-2
7.2	Procedure.....	7-3
7.3	Sulfur Measurement Kits.....	7-4
	Sulfur Measurement Kit for ConFlo III	7-4
	Sulfur Measurement Kit for Flash Elemental Analyzer.....	7-4

Finnigan ConFlo III

7.4	Preparing the System for a Sulfur Measurement	7-6
	Installing the Teflon Tubing	7-6
	Installing the Self-Adhesive Heating Foil	7-7
	Reactor Filling.....	7-8
7.5	Before Starting a Sulfur Measurement.....	7-9
	For EA 1108, EA 1110 and NA 2500.....	7-9
	For Flash Elemental Analyzer.....	7-10
	Stainless Steel GC Column.....	7-10
	Teflon GC Column	7-11
7.6	Create a Gas Configuration for a Sulfur Measurement.....	7-12
7.7	Zero Enrichment of Sulfur (Standard On/Off Test)	7-15
	Instrument.....	7-15
	Time Events.....	7-16
	Evaluation.....	7-17
	Peak Detection.....	7-18
	Printout	7-19
7.8	Linearity Test of Sulfur	7-22
Technical Information		8-1
8.1	Mechanics	8-2
8.2	Capillaries and Ferrules	8-4
8.3	Function Schematic.....	8-6
8.4	Connections for Compressed Air.....	8-7

Read This First

Welcome to the Thermo Electron, Finnigan ConFlo III Universal Interface!

This **Finnigan ConFlo III Operating Manual** describes how to setup and use the Finnigan ConFlo III Universal Interface.

This manual includes the following chapters:

Chapter 1: Prior to Measurement describes precaution measures and important steps to be performed in the beginning before a measurement is started.

Chapter 2: The ConFlo III Interface deals with the open split principle and treats the front and rear panel of ConFlo III respectively.

Chapter 3: ConFlo III Interface and Elemental Analyzer describes how to install an autosampler, how to operate ConFlo III manually and how to get started with ConFlo III plus Elemental Analyzer.

Chapter 4: Nitrogen Measurement describes zero enrichment (standard on/off test) and linearity test of nitrogen.

Note. This chapter treats the principles of blank measurement, reference measurement and amount percent determination. As they can be generalized for nitrogen, carbon and sulfur, the chapter already contains the remarks valid for any of the three elements.

Chapter 5: Carbon Measurement describes zero enrichment (standard on/off test) and linearity test of carbon.

Chapter 6: Dual Measurement describes the procedure and how to perform a jump calibration.

Chapter 7: Sulfur Measurement describes the sulfur measurement kit and how to prepare the system for a sulfur measurement. Besides, this chapter treats zero enrichment (standard on/off test) and linearity test of sulfur.

Chapter 8: Technical Information describes mechanics, capillaries, ferrules, the function schematic and connections for compressed air.

Changes to the Manual and Online Help

To suggest changes to this manual or the online Help, please send your comments to:

Thermo Electron (Bremen) GmbH
Finnigan Advanced Mass Spectrometry
Product Marketing
Barkhausenstr. 2

D-28197 Bremen

Germany

e-mail: documentation@thermo-bremen.com

You are encouraged to report errors or omissions in the text or index.

Thank you.

Abbreviations

The following abbreviations are used in this and other manuals and in the online Help.

A	ampere
ac	alternating current
ADC	analog-to-digital converter
AP	acquisition processor
APCI	atmospheric pressure chemical ionization
API	atmospheric pressure ionization
ASCII	American Standard Code for Information Interchange
b	bit
B	byte (8 b)
baud rate	data transmission speed in events per second
°C	degrees Celsius
CD	compact disc
CD-ROM	compact disc read-only memory
cfm	cubic feet per minute
CI	chemical ionization
CIP	carriage and insurance paid to
cm	centimeter
cm ³	cubic centimeter
CPU	central processing unit (of a computer)
CRC	cyclic redundancy check
CRM	consecutive reaction monitoring
<Ctrl>	control key on the terminal keyboard
<i>d</i>	depth
Da	dalton
DAC	digital-to-analog converter
dc	direct current
DDS	direct digital synthesizer
DEP™	direct exposure probe
DS	data system
DSP	digital signal processor

EI	electron ionization
EMBL	European Molecular Biology Laboratory
<Enter>	enter key on the terminal keyboard
ESD	electrostatic discharge
ESI	electrospray ionization
eV	electron volt
f	femto (10^{-15})
°F	degrees Fahrenheit
.fasta file	extension of a SEQUEST search database file
FOB	free on board
ft	foot
FTP	file transfer protocol
g	gram
G	giga (10^9)
GC	gas chromatograph; gas chromatography
GC/MS	gas chromatograph / mass spectrometer
GND	electrical ground
GPIB	general-purpose interface bus
GUI	graphical user interface
h	hour
<i>h</i>	height
HPLC	high-performance liquid chromatograph
HV	high voltage
Hz	hertz (cycles per second)
ICIS™	Interactive Chemical Information System
ICL™	Instrument Control Language™
ID	inside diameter
IEC	International Electrotechnical Commission
IEEE	Institute of Electrical and Electronics Engineers
in.	inch
I/O	input/output
k	kilo (10^3 , 1000)
K	kilo (2^{10} , 1024)
KEGG	Kyoto Encyclopedia of Genes and Genomes
kg	kilogram

<i>l</i>	length
L	liter
LAN	local area network
lb	pound
LC	liquid chromatograph; liquid chromatography
LC/MS	liquid chromatograph / mass spectrometer
LED	light-emitting diode
μ	micro (10^{-6})
m	meter
m	milli (10^{-3})
M	mega (10^6)
M+	molecular ion
MB	Megabyte (1048576 bytes)
MH+	protonated molecular ion
min	minute
mL	milliliter
mm	millimeter
MS	mass spectrometer; mass spectrometry
MS	MS ⁿ power: where n = 1
MS/MS	MS ⁿ power: where n = 2
MS ⁿ	MS ⁿ power: where n = 1 through 10
<i>m/z</i>	mass-to-charge ratio
n	nano (10^{-9})
NCBI	National Center for Biotechnology Information (USA)
NIST	National Institute of Standards and Technology (USA)
OD	outside diameter
Ω	ohm
p	pico (10^{-12})
Pa	pascal
PCB	printed circuit board
PID	proportional / integral / differential
P/N	part number
P/P	peak-to-peak voltage

ppm	parts per million
psig	pounds per square inch, gauge
RAM	random access memory
RF	radio frequency
RMS	root mean square
ROM	read-only memory
RS-232	industry standard for serial communications
s	second
SIM	selected ion monitoring
solids probe	direct insertion probe
SRM	selected reaction monitoring
SSQ [®]	single stage quadrupole
TCP/IP	transmission control protocol / Internet protocol
TIC	total ion current
Torr	torr
TSQ [®]	triple stage quadrupole
u	atomic mass unit
URL	uniform resource locator
V	volt
V ac	volts alternating current
V dc	volts direct current
vol	volume
w	width
W	watt
WWW	World Wide Web

Note. Exponents are written as superscripts. In the corresponding online Help, exponents are sometimes written with a caret (^) or with *e* notation because of design constraints in the online Help. For example:

MSⁿ (in this manual) Ms^{^n} (in the online Help)

10⁵ (in this manual) 10^{^5} (in the online Help)

Typographical Conventions

Typographical conventions have been established for Thermo Electron San Jose manuals for the following:

- Data input
- Boxed information
- Topic headings

Data Input

Throughout this manual, the following conventions indicate data input and output via the computer:

- Messages displayed on the screen are represented by capitalizing the initial letter of each word and by italicizing each word.
- Input that you enter by keyboard is represented in **bold face letters**. (Titles of topics, chapters, and manuals also appear in bold face letters.)
- For brevity, expressions such as “choose **File > Directories**” are used rather than “pull down the File menu and choose Directories.”
- Any command enclosed in angle brackets < > represents a single keystroke. For example, “press <F1>” means press the key labeled *F1*.
- Any command that requires pressing two or more keys simultaneously is shown with a plus sign connecting the keys. For example, “press <Shift> + <F1>” means press and hold the <Shift> key and then press the <F1> key.
- Any button that you click on the screen is represented in bold face letters and a different font. For example, “click on **Close**”.

Boxed Information

Information that is important, but not part of the main flow of text, is displayed in a box such as the one below.

Note. Boxes such as this are used to display information.

Boxed information can be of the following types:

- **Note** – information that can affect the quality of your data. In addition, notes often contain information that you might need if you are having trouble.
- **Caution** – information necessary to protect your instrument from damage.
- **Warning** – hazards to human beings. Each Warning is accompanied by a Warning symbol.

Topic Headings

The following headings are used to show the organization of topics within a chapter:

Chapter 1

Chapter Name

1.2 Second Level Topics

Third Level Topics

Fourth Level Topics

Fifth Level Topics

Reply Cards

Thermo Electron San Jose manuals contain one or two reply cards. All manuals contain a Customer Registration / Reader Survey card and some contain a Change of Location card. These cards are located at the front of each manual.

The Customer Registration / Reader Survey card has two functions. First, when you return the card, you are placed on the Thermo Electron San Jose mailing list. As a member of this list, you receive application reports and technical reports in your area of interest, and you are notified of events of interest, such as user meetings. Second, it allows you to tell us what you like and do not like about the manual.

The Change of Location card allows us to track the whereabouts of the instrument. Fill out and return the card if you move the instrument to another site within your company or if you sell the instrument. Occasionally, we need to notify owners of our products about safety or other issues.

Chapter 1

Prior to Measurement

1.1 Precautions



Warning. Ensure that ConFlo III is not exposed to direct sunlight!



Warning. Keep the system away from heat, e.g. radiators, to avoid damage to the internal circuits and the external surface!



Warning. Avoid using aerosol sprays near the system as this could cause sudden ignition of the spray!



Warning. Do not clean ConFlo III using paint thinner, alcohol or other organic solvents!



Warning. Do not put any objects, especially liquids, upon ConFlo III!



Warning. Place the system on a flat and solid surface that can carry at least 100 kg!

1.2 About this Manual

This manual intends to be a user guide for working with the combined system of Elemental Analyzer (EA), ConFlo III and Isotope Ratio Mass Spectrometer (IRMS). As the system consists of three separate instruments with the Elemental Analyzer and IRMS having manuals of their own, this manual will mark on specific information about the ConFlo III system. It starts with an overview about Elemental Analyzers of different types followed by basic information about ConFlo III. Finally, isotope ratio determinations of C, N, S, H and O using ConFlo III are explained in turn. It is assumed that the user is not only familiar with clean operating procedures and sample preparation but also has already some working experience with IRMS and the software Isodat 2.0.

1.3 Introduction

The Thermo Electron Continuous Flow Interface (ConFlo III) provides an efficient and flexible way to couple an Elemental Analyzer (EA) to a stable isotope ratio mass spectrometer (IRMS) on-line. This combination has become standard for determining the isotope ratios of carbon (C), nitrogen (N), sulfur (S), hydrogen (H) and oxygen (O) of combustible organic materials in bulk samples. ConFlo III and Elemental Analyzer form an integrated and optimized sample preparation inlet system, which can be connected to any current Thermo Electron IRMS equipped for on-line isotope ratio analysis. As a particularity of gas isotope mass spectrometers, any samples must be converted into simple gases prior to introducing them into the ion source. This is achieved by the Elemental Analyzer via oxidation and reduction furnaces (Elemental Analyzer) or by High Temperature Conversion (TC/EA). Passing through a gaschromatographic column the produced gases are separated and consecutively enter ConFlo III.

ConFlo III can be connected to any current Thermo Electron Elemental Analyzer and Thermo Electron IRMS equipped for Continuous Flow application.

Chapter 2

The ConFlo III Interface

2.1 Introduction - The Open Split Principle

The ConFlo III Interface provides the means for coupling the EA to the IRMS. The connection of an Elemental Analyzer with an isotope ratio mass spectrometer via an open split arrangement has become standard for on-line isotope ratio analysis. The IRMS works with a helium flow of approximately 0.3 ml/min, whereas the EA is operated with a helium flow of 80 to 120 ml/min. Reducing this gas flow is the central idea of the open split arrangement.

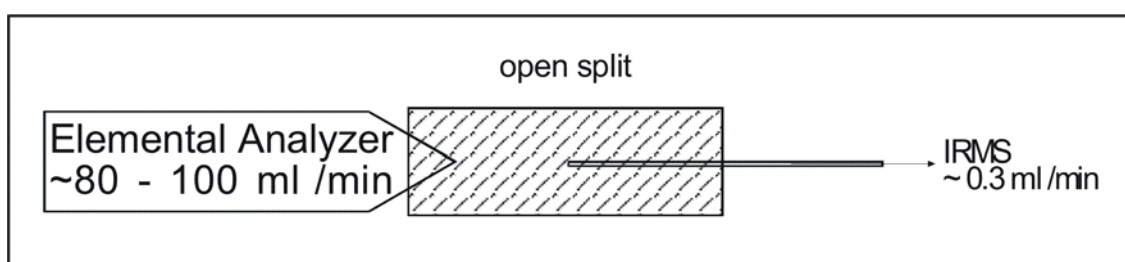


Figure 2-1. The Open Split Principle

The size of the IRMS capillary determines a gas flow of approximately 0.3 ml/min to the mass spectrometer (Figure 2-1). This IRMS capillary continuously "sniffs" the carrier gas (He) flowing through the cell. Hence, it may be referred to as the "sniffer" capillary. The carrier gas flushes the open split cell continuously with helium and passes the gaseous sample so transporting the produced compounds consecutively through the cell.

Procedure

The ConFlo III Interface forms an innovation in coupling an Elemental Analyzer to an IRMS in the way that two separate open split cells replace the "tube in tube" sample gas inlet performed in the predecessor model ConFlo II.

Through the separation of the sample stream and the reference inlet, ConFlo III forms a dynamic and flexible way of an EA/IRMS connection. These two open split cells, in the following titled as "sample section" and "reference section", allow a higher precision of on-line determination as well as a temporal independent sample/reference gas introduction to the IRMS (Figure 2-2).

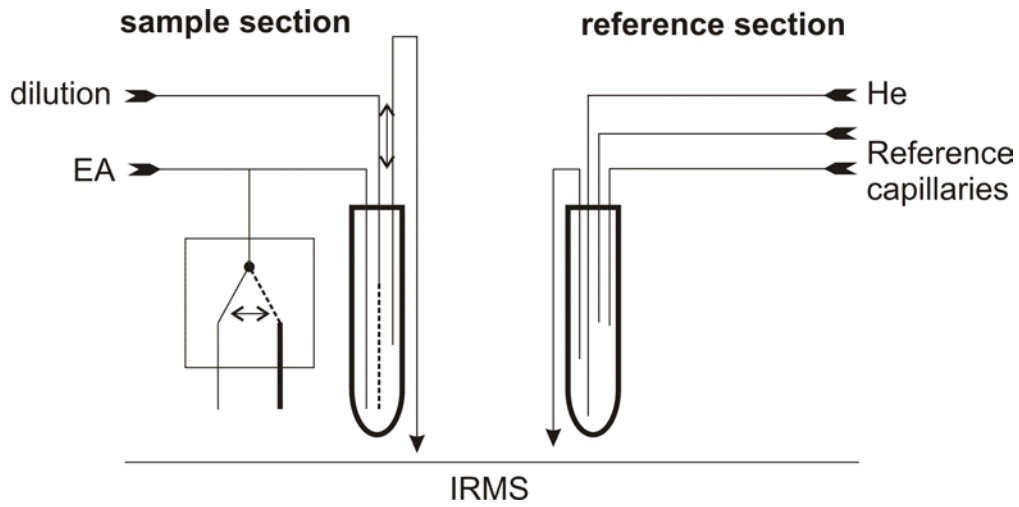


Figure 2-2. Sample Section and Reference Section

The Sample Section

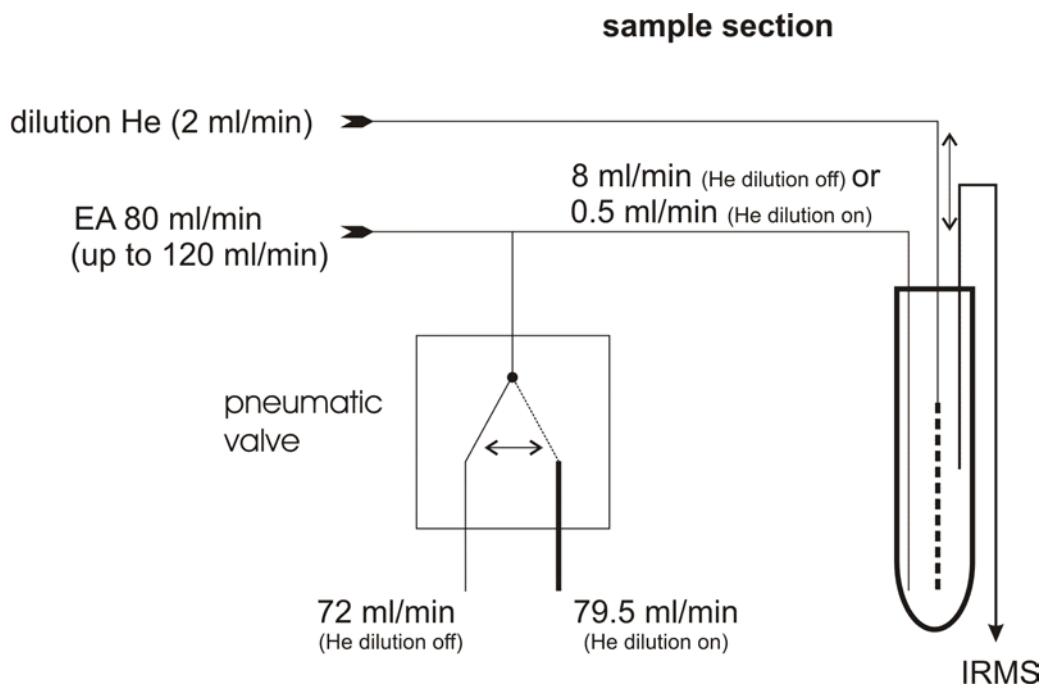


Figure 2-3. Sample Section

The open split arrangement of the sample section consists of three capillaries (Figure 2-3). The active, piston regulated dilution capillary, the fixed sample (delivering) capillary (which is located upstream the sniffer point) and the fixed "sniffer" (IRMS) capillary.

The effluent stream from the Elemental Analyzer is passed through a pneumatic valve before it enters the open split cell of the sample section. In case of dilution, the pneumatic valve is opened and the sample flow of the Elemental Analyzer is split in fixed relation (79.5 ml/min). Therefore, only 0.5 ml/min are delivered to the open split cell of the sample section. Simultaneously to the valve operation the piston with the dilution capillary moves down beyond the "sniffer" point. The He dilution flow can be varied manually between 5 and 50 ml/min resulting in an effect between 10 and 100-fold. The "sniffer" (IRMS) capillary with an inner diameter of 100 μm allows a flow of approximately 0.3 ml/min into the ion source. The dilution capillary is parked outside the "sniffed area", if dilution of the sample is not necessary.

The Reference Section

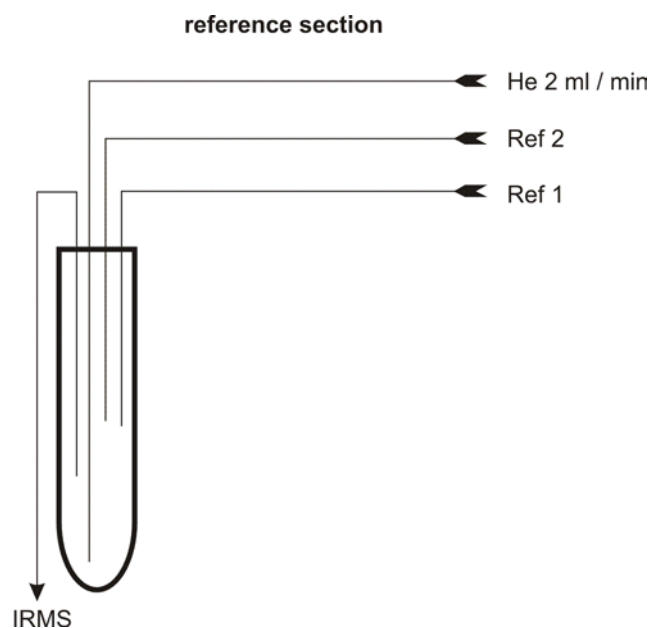


Figure 2-4. Reference Section

The open split cell of the reference section consists of four capillaries (Figure 2-4). The He carrier gas capillary as well as the IRMS capillary are fixed while a permanent stream of helium prevents a dry out of the IRMS capillary. In addition, two piston operated reference capillaries are located in the open split cell. Each reference capillary carries a continuous flow of gas

e.g. nitrogen and carbon dioxide. When a standard is required during an acquisition, the piston with the appropriate reference capillary moves down beyond the "sniffing" position and the standard is sniffed via the IRMS capillary to the mass spectrometer. Because the standard gas is injected directly into the open split cell, the mass spectrometer response is almost instantaneous, delayed only by the time the gas needs to pass through the capillary.

2.2 Front and Rear Panel of ConFlo III

Front Panel

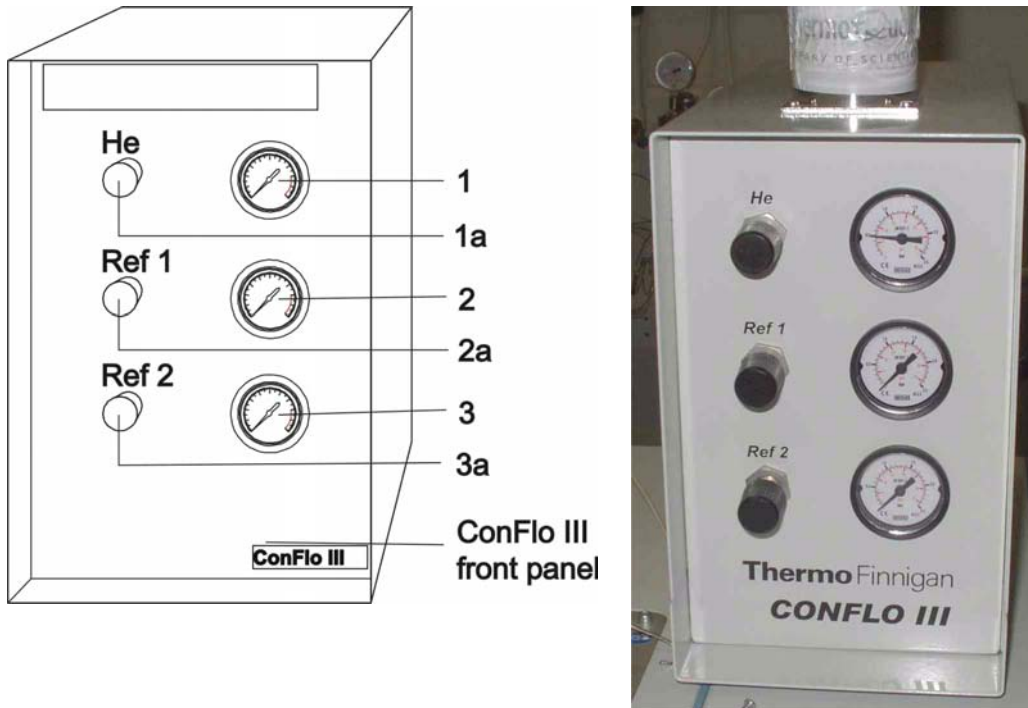


Figure 2-5. Front Panel of ConFlo III

- 1) Helium pressure gauge
- 1a) Helium pressure regulator
- 2) Reference gas 1 pressure gauge
- 2a) Reference gas 1 pressure regulator
- 3) Reference gas 2 pressure gauge
- 3a) Reference gas 2 pressure regulator

Side Views

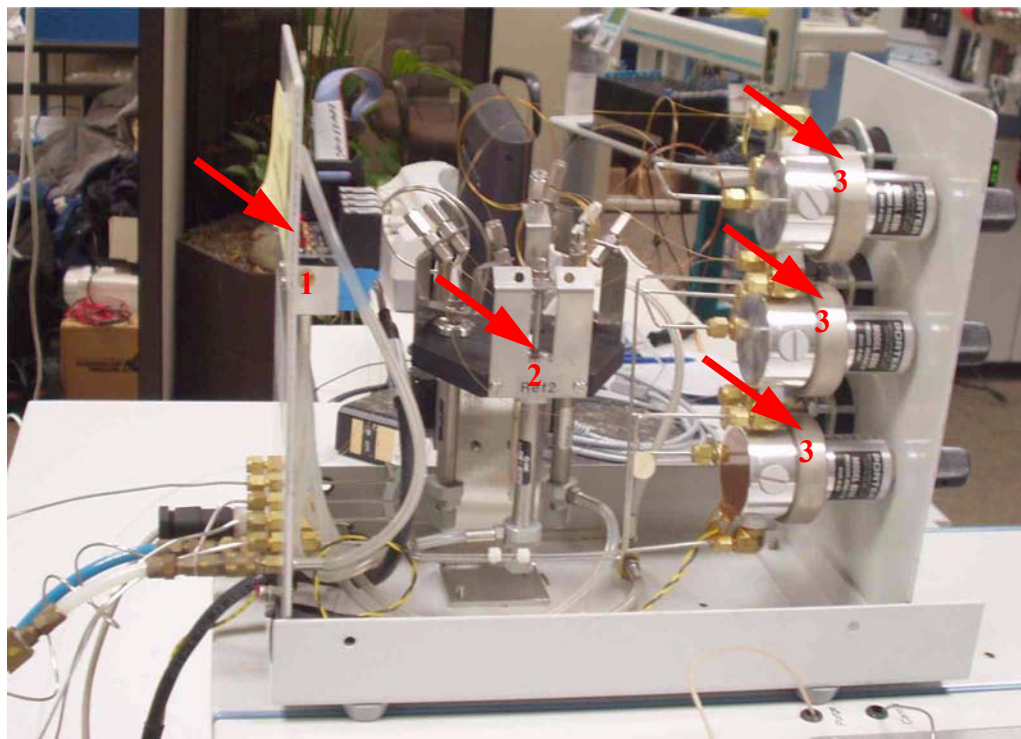


Figure 2-6. ConFlo III - Left Side View

- 1 manifold card
- 2 combi inlet with piston-operated actuators
- 3 pressure gauge

For technical information see also **Chapter 8: Technical Information.**

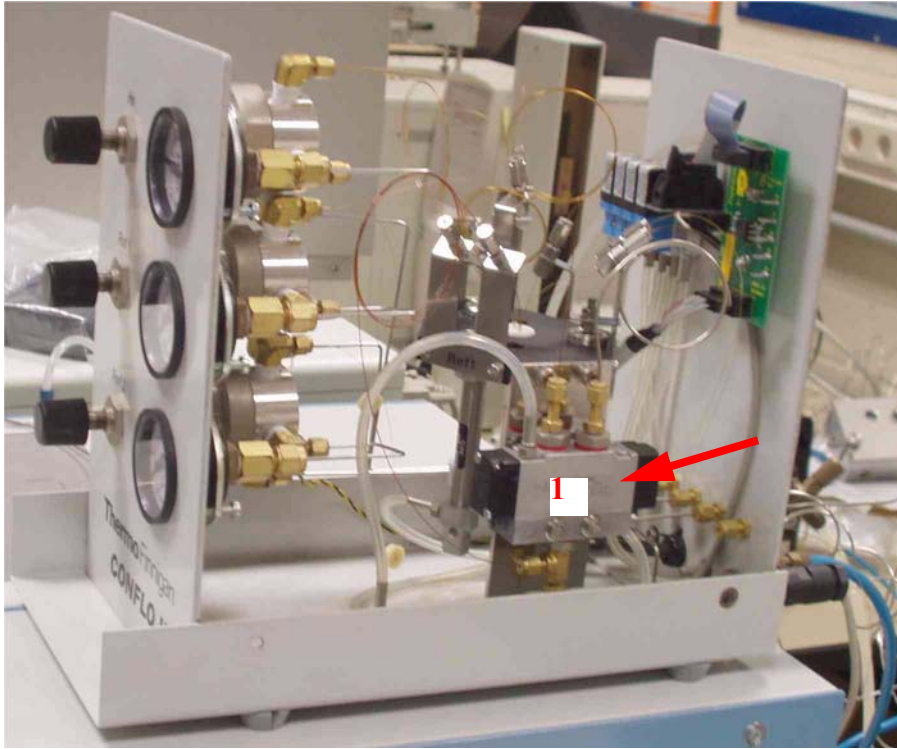


Figure 2-7. ConFlo III - Right Side View

- 1 two port valve (prediluter); see also Figure 8-1.

Rear Panel and Gas Connections

External gas connections are located at the rear panel of ConFlo III (Figure 2-8).

1. Fused silica port for the IRMS capillary
2. Remote Elemental Analyzer start cable
3. Compressed air outlet (time controlled)
4. Elemental Analyzer inlet
5. Reference gas 1 inlet
6. Reference gas 2 inlet
7. Helium inlet (carrier gas)
8. Compressed air inlet (permanent)
9. IRMS connection cable

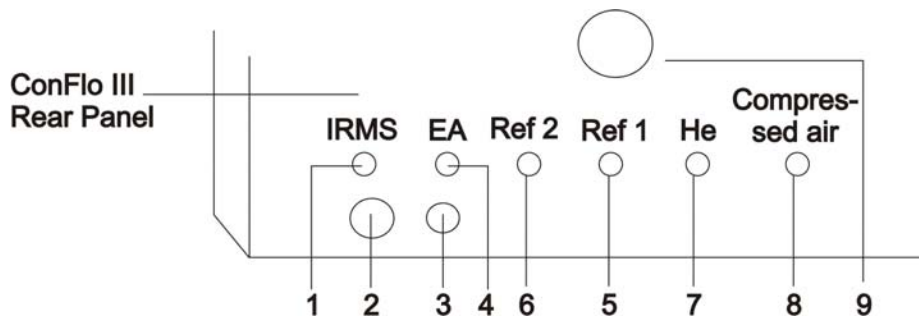


Figure 2-8. Rear Panel



Warning. For detailed information, refer to the gas flow diagram (Figure 2-9). Wrong handling of the toxic gases CO and SO₂ is perilous! Inform yourself how to handle these gases by reading the directions for use and asking your local gas tank seller. Thermo Electron does not take any responsibility if the standard gas (CO) or other gases, e.g. SO₂, are used incorrectly.

Gas Flow Diagram

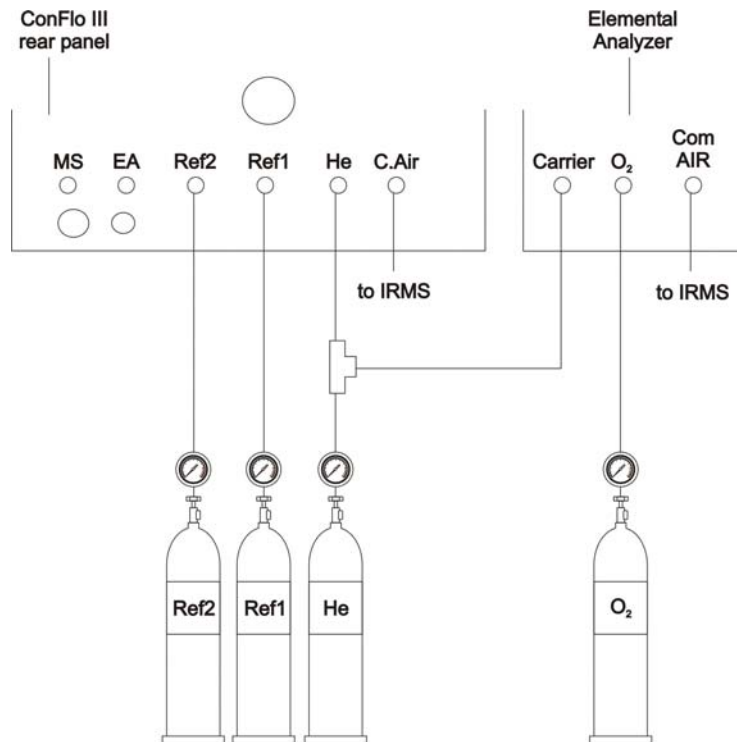


Figure 2-9. Gas Flow Diagram

Spare Parts of Reference and Sample Sections

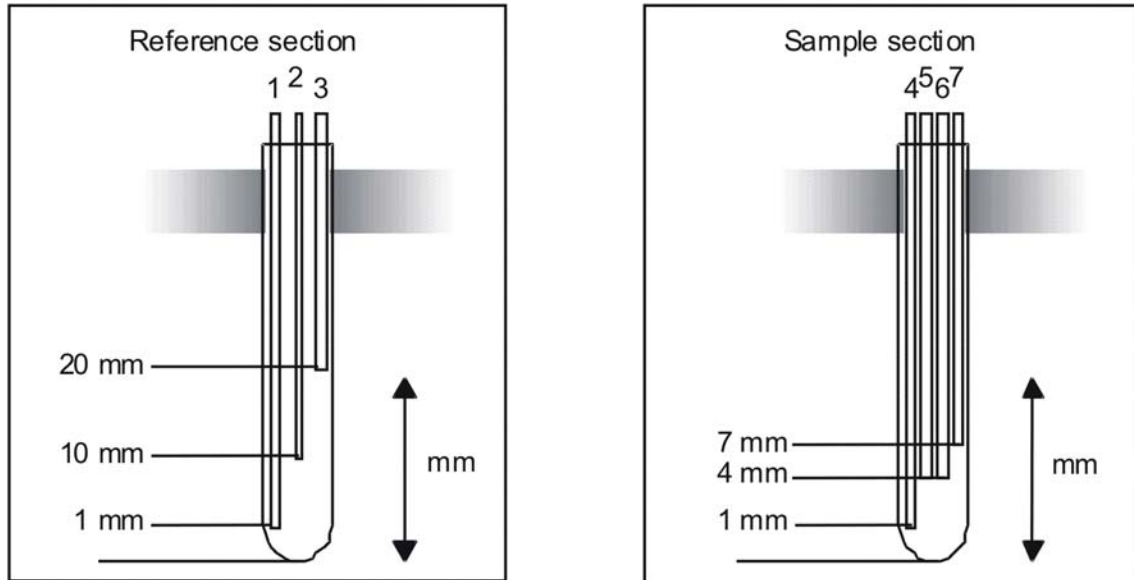


Figure 2-10. Reference Section and Sample Section

Table 2-1. Spare Parts of Reference and Sample Sections - Part 1

Capillary	Length	Inner diameter	Part No.
1 Fused silica from EA (T-piece)	1000 mm	0.32 mm	100 4640
2 Fused silica helium dilution	2000 mm (or 1000 mm)	0.18 mm	054 2910
3 Fused silica to IRMS	2000 mm	0.10 mm	104 8990
4 Fused silica helium purge	250 mm (or 750 mm)	0.10 mm	104 8990
5 Fused silica Ref. 1	310 mm (or 2000 mm)	0.05 mm	054 3380
Vent	200 mm	0.075 mm	104 5480
6 Fused silica Ref. 2	340 mm (or 2000 mm)	0.05 mm	054 3380
Vent	200 mm	0.075 mm	104 5480
7 Fused silica to IRMS	200 mm	0.10 mm	104 8990
	1500 mm	0.05 mm	054 3380

Table 2-2. Spare Parts of Reference and Sample Sections - Part 2

	Capillary	Inner diameter	Ferrule	Part No.
1	Fused silica	0.32 mm	single hole	056 6390
2	Fused silica	0.18 mm	single hole	100 4850
3	Fused silica	0.10 mm	single hole	100 4850
4	Fused silica	0.05 mm	single hole	100 4850

Manometer

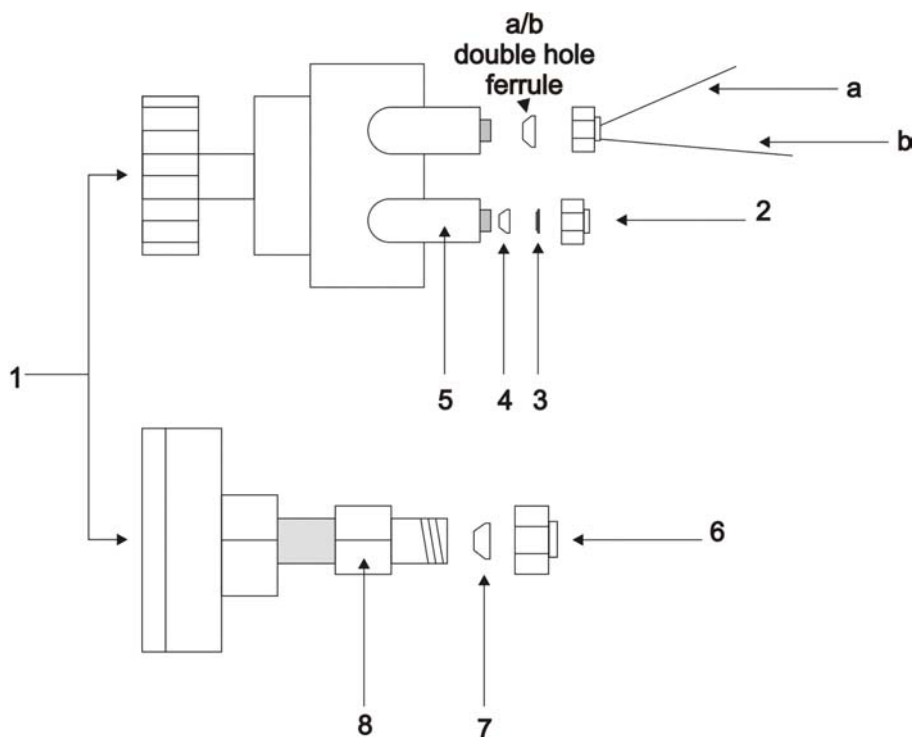


Figure 2-11. Manometer

Table 2-3. Manometer Parts

Pos.	Designation	Part No.
a)	He dilution	a / b double hole ferrule: 106 0170
b)	He purge	
a)	Reference gas	a / b double hole ferrule: 100 4780
b)	Reference purge	
1	Manometer 0 - 2.5 bar	052 43 91
2	Cap nut 1/16"	052 3610
3	Back ferrule 1/16"	052 3210
4	Front ferrule 1/16"	052 3440
5	Elbow	049 3190
6	Cap nut 1/8"	052 1500
7	Ferrule; Teflon 1/8"	067 4790
8	Coupling 1/8"	074 3360

Chapter 3

ConFlo III Interface and Elemental Analyzer

3.1 ConFlo III Software in Old Isodat

The ConFlo III application software allows fully automated determination of the isotope ratios of carbon, nitrogen, oxygen, hydrogen and sulfur from combustible organic matter.

Note. The ConFlo III software is used within the old Isodat software. Therefore, it is necessary to have the old Isodat software properly installed.

The ConFlo III software is installed as follows:

- Stop computer from *Isodat* mode by holding the *Alt* key while typing *Quit*.
- Press *Ctrl* and *4* simultaneously. The computer shows: c:\>ISOTST.
- Type: c:\>ISOTST>*install*.



Note. If the message **Not enough memory** appears:
Type: **BANK=ON**.
Press **ENTER**.

- Press *ENTER*. The <*Isodat*> *Installation Menu* appears:
 1. GEM Installation Menu
 2. ...
 3. ...
 6. GC/ConFlo III Installation Menu
 7. ...
 - Press **6**, then *ENTER*.
 - The *GC/ConFlo Installation Menu* appears:
 1. ...
 2. ConFlo III Installation
 3. ...
 - Press **2**, then *ENTER* and continue the procedure.
 - After finishing boot the system by pressing *Ctrl+Alt+Del* simultaneously.

3.2 ConFlo III Configuration in Isodat 2.0



Configurator

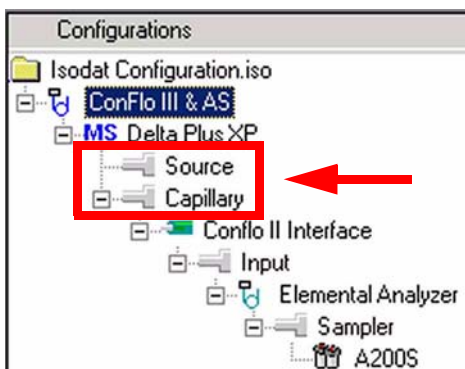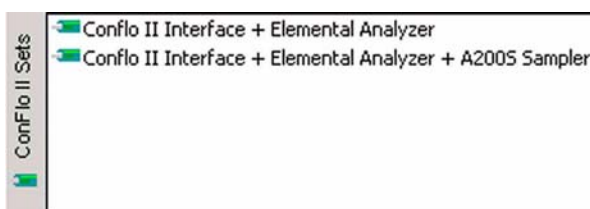

- Before operating, a **Configuration** containing the ConFlo II/III Interface needs to be created in the **Configurator**.
 - Press the **Add Configuration** button to add a new Configuration.
 - Give it a name, e.g. "ConFlo III & AS".
 - Select one of the ConFlo II sets (also for ConFlo III) and drag it either to the **Source** port or to the **Capillary** port in the new Configuration.
 - In case of a ConFlo Interface and an autosampler with a RS 232 interface to the computer select **ConFlo Interface + Elemental Analyzer + A200S Sampler**.
 - Open the complete tree structure of the ConFlo Interface to check for the hardware attached (Elemental Analyzer, autosampler).
 - In case of a ConFlo II/III and a liquid autosampler with a RS 232 interface to the computer, the COM Port settings can be checked in the **Advanced Mode**.
- See **Installation of the A200S and GC PAL Autosampler** on page 3-4.
- Close the Configurator window. All settings will be saved automatically.


3.3 Installation of the A200S and GC PAL Autosampler

A200S and GC PAL autosampler require operating signals from the software via RS 232 cable (COM Port). Special scripts for using these autosamplers together with the software Isodat 2.0 are installed automatically by the selection of a ConFlo III set with liquid autosampler in the Isodat 2.0 *Configurator*.

The COM Port settings for the autosampler can be viewed using the *Advanced Mode* (i.e. *Edit > Advanced Mode*):



- In the *Edit* Menu, activate the *Advanced Mode*.

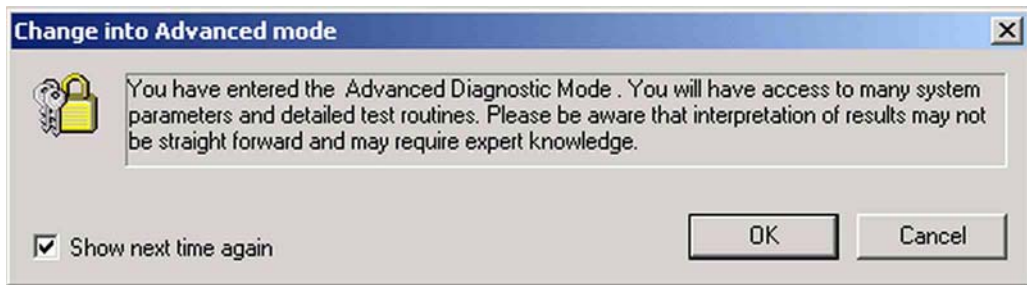


- Confirm by *OK*.

Note. The COM Port settings are essential for proper communication with the autosampler. Incorrect settings will lead to inactive liquid autosampler control.



- Press the *Devices* tab.

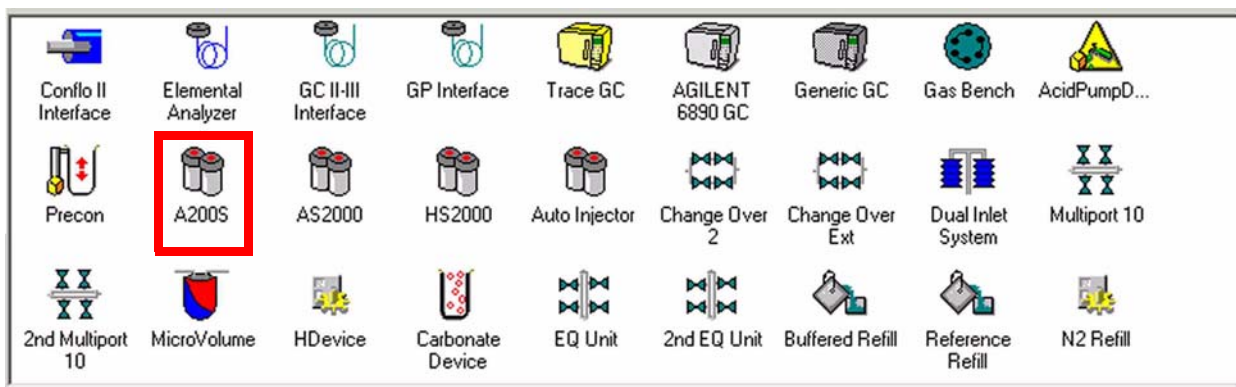

- The pool of available devices is shown.
- **Right-click** on your autosampler (e.g. A200S). For GC PAL autosampler also right-click A200S.
- Select **Properties**, if you only want to check the COM Port your autosampler is connected to or
- Select **Edit...** to edit the default communication parameters.







- If you selected **Properties** above, choose the COM Port your autosampler is connected to, e.g. COM 1.
- Confirm by **OK**.

Note. The **Properties** command is available also in **Normal Mode**, i.e. you don't need to be in **Advanced Mode** to only check the COM Port your autosampler is connected to.


The **Edit...** command, however, is available only in **Advanced Mode**.

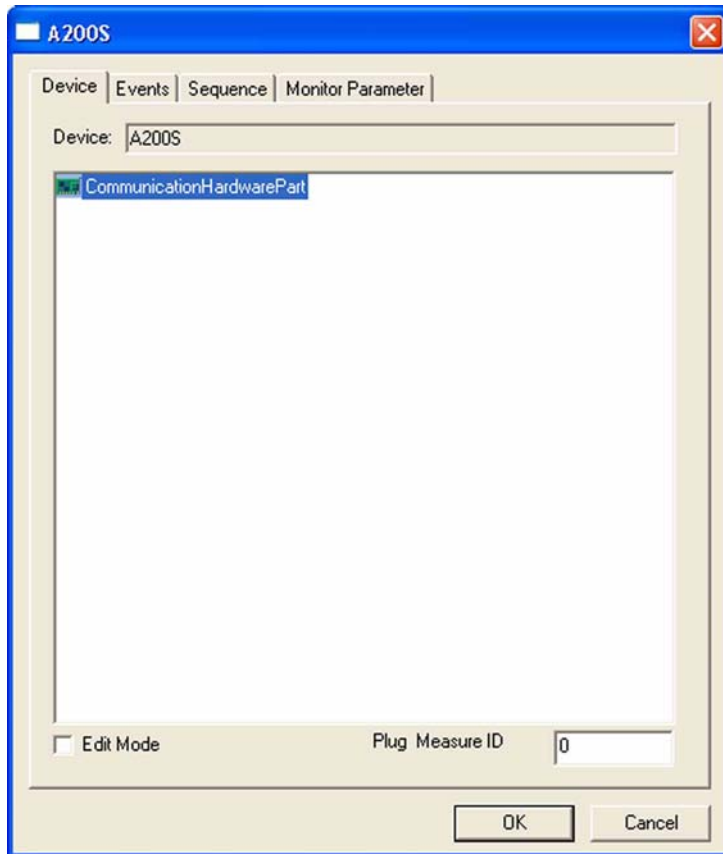

- If you selected *Edit...* above, double-click on *Communication Hardware Part*. Alternatively, right-click on it and select *Properties*.

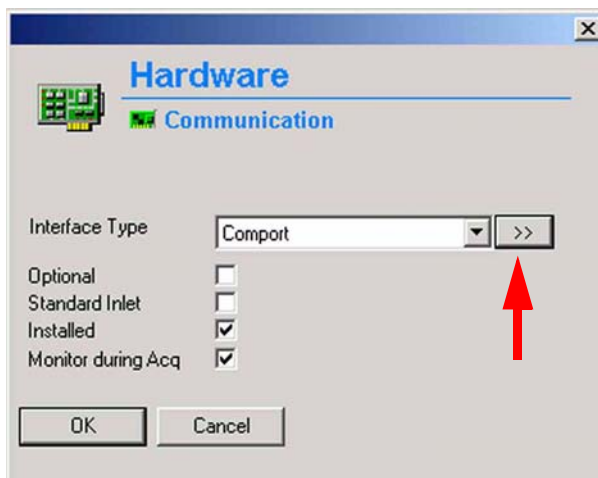

- The default settings are shown above. Usually, they need not to be edited.
- Choose your interface type (*Comport*).
- To specify its properties, press the >> button.

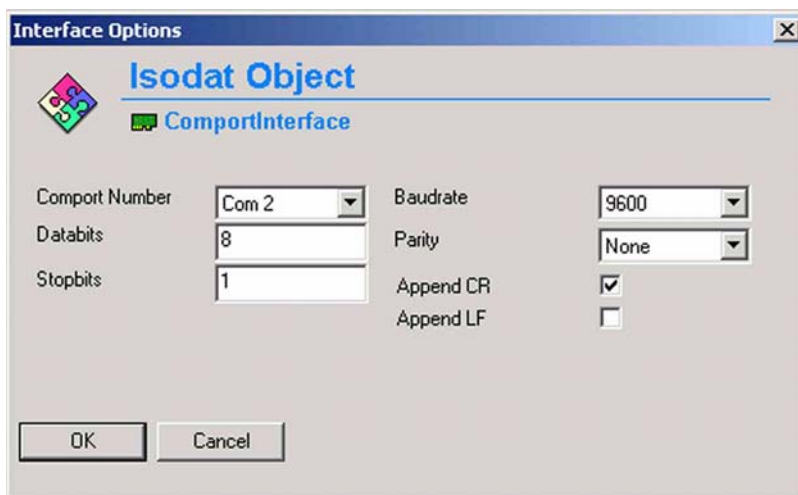

- The default settings are shown above (usually kept default).
- Check, whether the COM Port corresponds to the COM Port your autosampler is connected to. Confirm by **OK**.

Note. The COM Port settings are essential for proper communication. Incorrect settings will lead to inactive hardware control.


3.4 Manual Operation of ConFlo III



- Open the *Instrument Control* module.



- Select a *Configuration* that contains the ConFlo III interface.

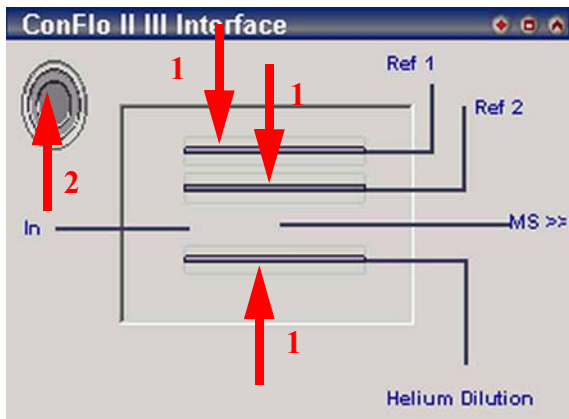

The *ConFlo II III Interface window* allows:
reference gas inlet and/or dilution (1)
autosampler/Elemental Analyzer start signal (2)

Capillaries can be moved/activated by mouse clicks.

- Open *Reference 1* by a *click* on *Ref 1* capillary. If CO₂ is connected to *Reference 2*, click *Ref 2* capillary.


Thus, the blue capillary symbol elongates and CO₂ enters the system.

Figure 3-1. ConFlo II III Interface Window

Comment: If the ConFlo II III Interface Window is not Visible



- *Right-click* somewhere into the *Accessories* toolbar.
- Click on the *Administrative Panels...* button.

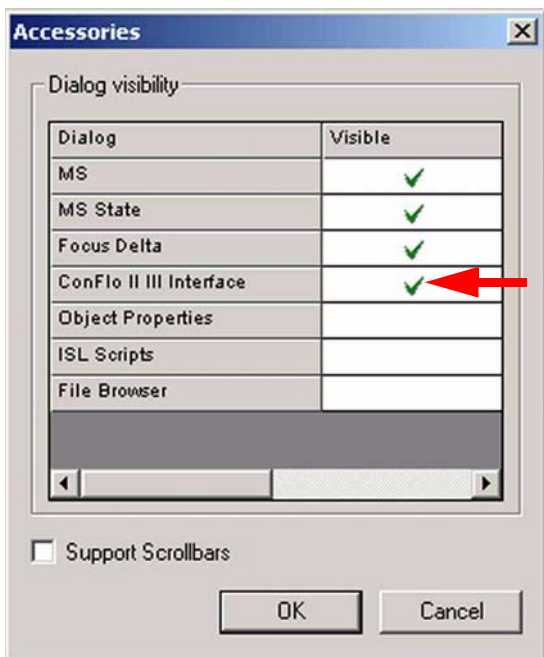

The *Accessories* window allows to mark dialogs (i.e. windows) by ✓. If you mark a dialog, the respective window will be shown at the *Accessories* toolbar.


- Mark *ConFlo II III Interface*.
- Confirm by **OK**.

Thus, the ConFlo II III Interface window will be shown at the *Accessories* toolbar.

Figure 3-2. Visible Windows at the Accessories Toolbar

Comment: If only the Title Bar of the ConFlo II III Interface Window is Visible



- Click on the *Unshrink* button  at the right corner of the Title bar.
Thus, ConFlo II III Interface window will appear.

3.5 How to Get Started with ConFlo III and Elemental Analyzer

1. Connect the Elemental Analyzer to ConFlo III for the transfer of the start signal.
2. In case you have Teflon tubing, replace it with stainless steel at your Elemental Analyzer, because atmospheric N₂ passes Teflon and interferes the quality of the determination.
3. Connect the reducing valves to the gas supplies (He, O₂, N₂, CO₂).
4. Connect the gas tank of O₂ and He to the Elemental Analyzer.
5. Connect He, N₂ (as Ref 2) and CO₂ (as Ref 1) to ConFlo III.
6. Open the main valves on the gas tanks and adjust the pressure.
7. Leak-check all connections and valves outside ConFlo III.
8. Make sure that the Elemental Analyzer is set up (i.e. oxidation and reduction reactors are packed and installed, water trap is installed, autosampler is fixed etc. Also refer to the Elemental Analyzer's manual).
9. Connect the Elemental Analyzer to the computer, if the Elemental Analyzer is run by separate software, i.e. Eager software.
10. Connect the Elemental Analyzer to the 220 V main power supply.
11. Connect compressed air to ConFlo III.
12. Connect compressed air to Elemental Analyzer if necessary (depending on the autosampler).
13. Connect ConFlo III to the IRMS (rear panel).
14. Ensure that the IRMS connection cable is connected to the driver board of the computer.
15. Make sure that the ConFlo III configuration has been installed during Isodat 2.0 setup.
See **ConFlo III Configuration in Isodat 2.0** on page 3-3.
16. By using the ConFlo III device window in Isodat 2.0, make sure that the respective pistons of ConFlo III are movable (click on them using your mouse) and watch them moving in the combi inlet of ConFlo III.
17. Connect the fused silica capillaries of ConFlo III to the needle valve of the IRMS.
18. Switch on the Elemental Analyzer.
19. Adjust the gas pressures of the Elemental Analyzer and ConFlo III.
Take the following settings as a guideline:

Table 3-1. Adjustment of Gas Pressures

He	Elemental Analyzer	100 kPa
O ₂	Elemental Analyzer	120 kPa
Purge with	He	110 - 200 ml/min depending on your autosampler:
		<ul style="list-style-type: none"> • AS 200 and AS 128: 110 ml/min • MAS 200: 200 ml/min
He	ConFlo III	0.8 bar
N ₂	ConFlo III	1.3 bar
CO ₂	ConFlo III	0.9 bar

Caution. To prevent atmospheric air from entering the open split cell, never set the He flow on ConFlo III to less than 0.5 bar.

20. Switch off the ion source.
21. Inject the He dilution capillary to ensure only He entering the source.
22. Open the needle valve slowly.
23. Make sure the ion source high vacuum is between $3 * 10^{-6}$ mbar and $1.5 * 10^{-6}$ mbar.
24. Switch the ion source on. The expected background after approximately 30 min should be as described below:

Table 3-2. Expected Background After Approximately 30 min

mass	measured in cup for mass	resistor [Ω]	expected background
28	28	$3 * 10^8$	< 20 mV
29	29	$3 * 10^{10}$	< 20 mV
18	29	$3 * 10^{10}$	< 10 mV
40	29	$3 * 10^{10}$	< 10 mV

Note. Background values may vary depending on sensitivity and focus settings. They are given here as a general guideline.

25. Leak-check the needle valve.

Note. Leak check can be performed using a triggering stream of Ar while monitoring mass 40 on the IRMS.

- 26. Switch the needle valve heater **ON**.
- 27. Switch the ion source heater **ON**.
- 28. Leak-check all connections of the Elemental Analyzer following the procedure described in the respective Elemental Analyzer manual.
- 29. Heat up the following parts of the Elemental Analyzer in three steps.

Note. The Elemental Analyzer should not be connected to ConFlo III yet!

Table 3-3. Heating Up Elemental Analyzer Parts in Three Steps

	step 1	step 2	step 3
oxidation furnace	400 °C	750 °C	1020 °C
reduction furnace	400 °C	500 °C	650 °C
column	40 °C	75 °C	120 °C

- 30. Leak-check after each step and wait for two hours after reaching the final temperature.
- 31. Keep the system overnight in the last status to remove organic material and water and for column conditioning.
- 32. Also leave ConFlo III under above conditions overnight to remove all organic material from the fused silica capillaries, connection lines, valves, manometers etc.
- 33. Cool down the GC column to 40 °C for measurements.
- 34. Before connecting the Elemental Analyzer to ConFlo III, measure the He flow at the end of the stainless steel capillary.
Adjust the pressure at the Elemental Analyzer to set the flow to 80 ml/min - 100 ml/min depending on sample and application to be performed.
- 35. Connect the Elemental Analyzer to ConFlo III.
- 36. Make sure that no standard gas is injected, He dilution is **off** and all other valves are closed.

The expected background is:

Table 3-4. Expected Background

mass	measured in cup for mass	resistor [Ω]	expected background
28	28	$3 * 10^8$	< 20 mV
29	29	$3 * 10^{10}$	< 20 mV
18	29	$3 * 10^{10}$	< 2 mV
40	29	$3 * 10^{10}$	< 10 mV

Note. Background values may vary depending on sensitivity and focus settings. They are given here as a general guideline.

Note. The above ion intensity depends on the sensitivity of the IRMS. If the ion intensity is higher than expected and no leak is present, the background must decrease after two or three days and stabilize.

Chapter 4

Nitrogen Measurement

4.1 Zero Enrichment of Nitrogen (Standard On/Off Test)

We assume that the user already has working experience with the ConFlo III Interface and IRMS. It is recommended to perform a simple check in order to test the analytical condition of ConFlo III and IRMS before measuring any samples. The most important checks to test the analytical condition are *zero enrichment* and *linearity test*.

1. Only as a guideline use the following method from the File Browser's *Examples* folder (*N2_zero.met*):

Instrument

The screenshot shows the 'Instrument' configuration tab. The 'Reference Port' dropdown menu is set to 'Reference 2', indicated by a red arrow. Other visible settings include: Experiment: Continuous flow; Configuration: ConFlo; Gasconfiguration: N2; Acquisition Script: Acquisition.sct; Isotope MS: Integration Time: 0.200 [s]; Peak Center Cup: Cup 3; Peak Center Predelay (s): 20; Peak Center Postdelay (s): 10; Reference Device: Use Scripts (unchecked); Reference Port: Reference 2.

Figure 4-1. Zero Enrichment of Nitrogen - Instrument Tab

- Select the reference port your reference gas is connected to at the ConFlo interface (e.g. Reference 2).
- Adapt the reference gas port to the respective column of the Time Events list: e.g. if you choose Reference 2, the on-off entries must occur in the Reference 2-On column of the Time Events list (see Figure 4-2).

Time Events

Instrument	Time Events	Component Names	Evaluation@N2	Peak Detection@N2	Printout@N2	
Time [s]	Start Sampler	Elemental Anal. - On	Dilution - On	Reference 1 - On	Reference 2 - On	Switch Gas
20					●	
40						●
60					●	
80						●
100					●	
120						●
140					●	
160						●
180					●	
200						●
220					●	
240						●
260					●	
280						●
300					●	
320						●

Figure 4-2. Zero Enrichment of Nitrogen - Time Events Tab

- Recognize the eight on-off pulses shown above in the expected chromatogram (Figure 4-8).
- Note that the off-time of the e.g. fourth reference gas pulse is 160 s (Figure 4-4).

Acquisition Start	Immediately	Acquisition End Time [s]	350
-------------------	-------------	--------------------------	-----

Figure 4-3. Zero Enrichment of Nitrogen - Acquisition Start and Acquisition End Time

Evaluation

Instrument | Time Events | Evaluation@N2 | Peak Detection@N2 | Printout@N2

Evaluation Type: N2

Ref. Nr.:	Ref. Time:	Ref. Name:	d 15N/14N	vs.
1	160.00	N2_zero	0.000	Air-N2

Reference/Blank

Significant Peak Start [s]: 0.000 Significant Peak Stop [s]: 0.000

Amount Percent [%]: 0.000 Unit: mg

Figure 4-4. Zero Enrichment of Nitrogen - Evaluation Tab

Note. At **Ref. Time**, the **off**-value of an arbitrary reference gas peak according to the Time Events list must be typed (e.g. 160 s as the **off**-value of the forth reference gas peak, see Figure 4-2).

Peak Detection

Instrument | Time Events | Evaluation@N2 | **Peak Detection@N2** | Printout@N2

Perform Peak Detection Perform Background Detection Detection on Mass

Detection Parameter

Start Slope [mV/s]

End Slope [mV/s]

Peak Min Height [mV]

Peak Resolution [%]

Max Peak Width [s]

Perform Timeshift

Background Parameter

Background Type

History [s]

Auto Square Pulse Recognition / Timeshift Suppression

Enable Factor rArea / Pk Width / Pk Height

Advanced Parameter >>

Figure 4-5. Zero Enrichment of Nitrogen - Peak Detection Tab

<< Advanced Parameter

Peak Detection Parameter Sets

Nr.:	Start Detection [s]	Stop Detection [s]
1	-1.000	-1.000

Smoothing

Smooth Type Number Of Datapoints

Figure 4-6. Zero Enrichment of Nitrogen - Advanced Parameters in Peak Detection Tab

Note. A value of -1 denotes unlimited.

Printout

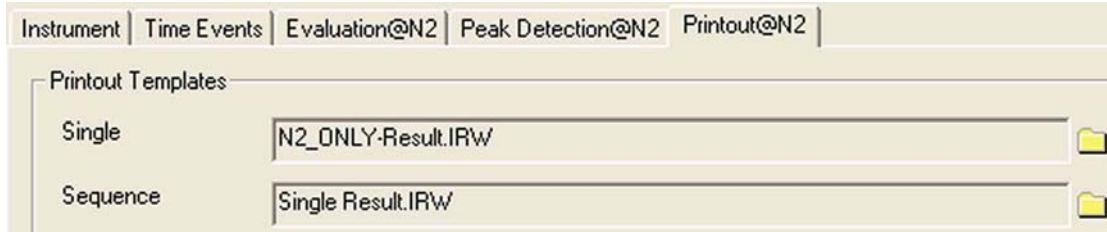


Figure 4-7. Zero Enrichment of Nitrogen - Printout Tab

2. At the corresponding pressure regulator of ConFlo III, set the ion intensity of mass 28 (i.e. $^{14}\text{N}^{14}\text{N}$) to 3 V - 4 V.
3. Create a new Sequence.
4. Press the **Start** button and confirm by **OK**.
5. Expected data after three or four measurements:

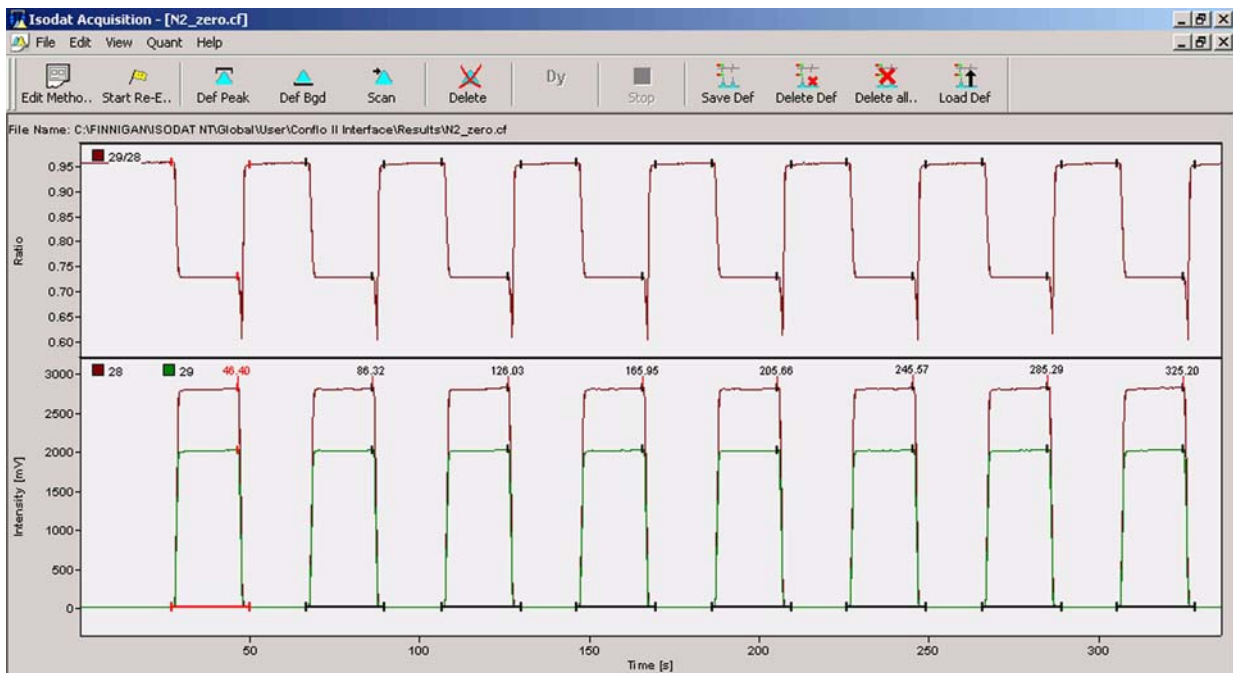


Figure 4-8. Zero Enrichment of Nitrogen - Chromatogram

Recognize the eight peaks shown above as the eight on-off pulses in the Time Events list (Figure 4-2).

N2		Error	Extended	Sequence Line						
Peak Nr.	Start [s]	Rt [s]	Width [s]	Ampl. 28 [mV]	Ampl. 29 [mV]	BGD 28 [mV]	BGD 29 [mV]	Area All [Vs]	d 15N/14N [per mil] vs. Air-N2	AT% 15N/14N [%]
1	27.0	46.4	23.2	2805	2016	17.0	12.0	52.610	-0.013	0.366467
2	66.9	86.3	23.0	2805	2017	17.5	12.4	52.656	0.004	0.366473
3*	106.6	126.0	23.4	2812	2022	17.6	12.5	52.599	0.000	0.366472
4	146.5	165.9	23.2	2810	2020	17.8	12.5	52.725	0.035	0.366485
5	186.2	205.7	23.4	2805	2016	17.7	12.5	52.696	0.036	0.366485
6	226.1	245.6	23.2	2815	2024	17.8	12.5	52.822	0.058	0.366493
7	265.8	285.3	23.4	2813	2023	17.8	12.6	52.720	0.037	0.366485
8	305.8	325.2	23.0	2810	2020	17.8	12.6	52.829	0.047	0.366489

Figure 4-9. Zero Enrichment of Nitrogen - Result Grid

Note. The standard deviation of $\delta^{15}\text{N}$ ($^{15}\text{N}^{14}\text{N}/^{14}\text{N}^{14}\text{N}$) should be less than 0.05 %.


To obtain the standard deviation of all eight peaks perform as follows:

d 15N/14N [per mil] vs. Air-N2
-0.013
0.004
0.000
0.035
0.036
0.058
0.037
0.047

- Click on the column header of the *d 15N/14N [per mil] vs. Air-N2* column. It will be highlighted.



- Right-click on the column header.
- Choose *Calculate*.

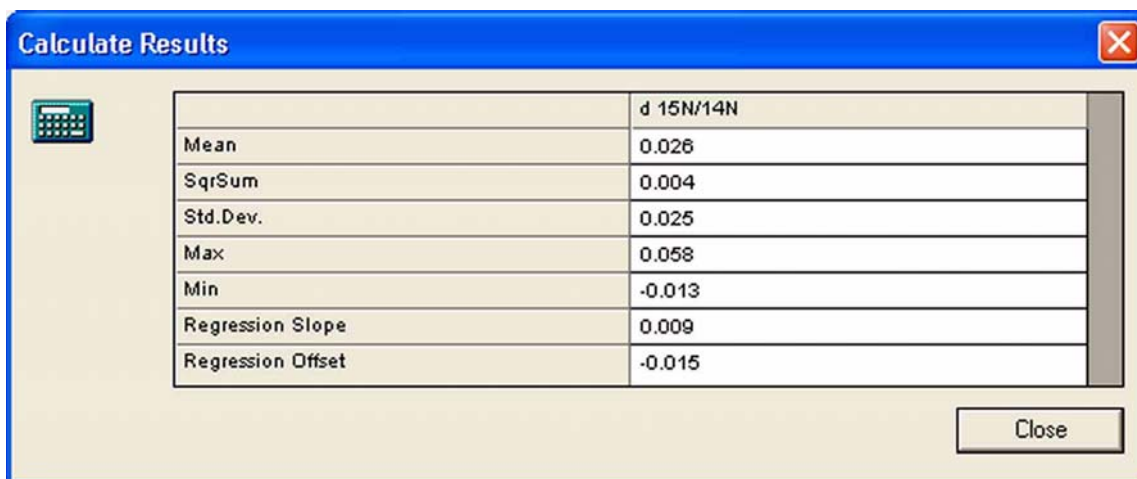

4.2 Linearity Test of Nitrogen

- Use the same method as defined for zero enrichment (see chapter **Zero Enrichment of Nitrogen (Standard On/Off Test)** on page 4-2).
- Start the acquisition as a single run.
- At each detection of a peak de- or increase the reference gas pressure at ConFlo III. Expected data after three or four measurements:

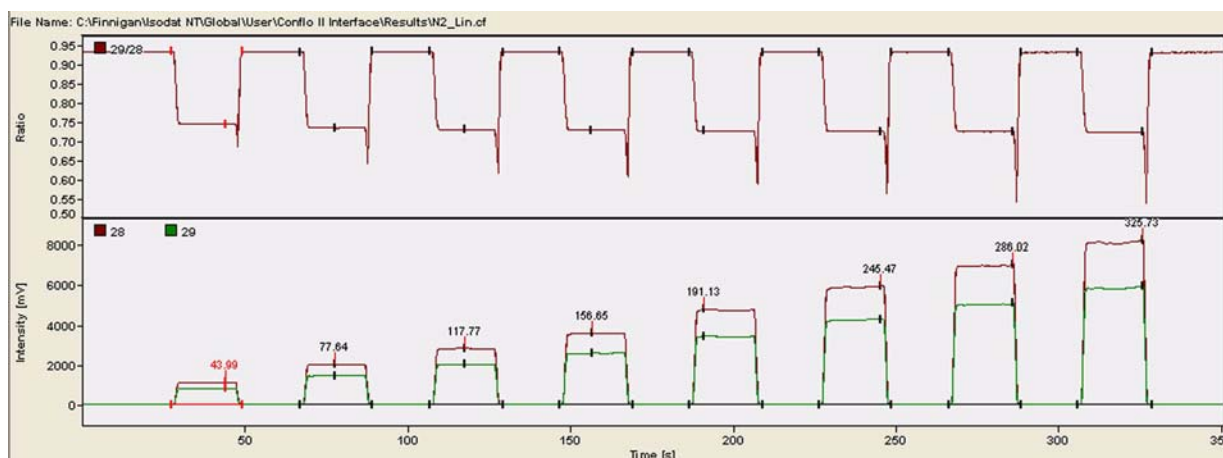


Figure 4-10. Linearity Test of Nitrogen - Chromatogram

N2	Error	Extended	Sequence Line							
Peak Nr.	Start [s]	Rt [s]	Width [s]	Ampl. 28 [mV]	Ampl. 29 [mV]	BGD 28 [mV]	BGD 29 [mV]	Area All [Vs]	d 15N/14N [per mil] vs. Air-N2	AT% 15N/14N [%]
1	27.5	44.0	21.9	1109	800	36.8	27.8	20.617	0.050	0.366490
2	67.2	77.6	22.2	2008	1449	36.7	27.7	37.623	0.101	0.366509
3*	106.9	117.8	22.4	2809	2026	36.8	27.7	52.679	0.000	0.366472
4	146.8	156.6	22.4	3577	2581	36.9	27.8	67.121	0.016	0.366478
5	186.5	191.1	22.6	4758	3433	37.0	27.9	89.271	0.046	0.366489
6	226.5	245.5	22.4	5932	4281	37.1	28.0	111.019	0.097	0.366507
7	266.4	286.0	22.4	7029	5072	37.3	28.1	131.249	0.113	0.366513
8	306.1	325.7	22.8	8194	5913	37.5	28.2	152.882	0.170	0.366534

Figure 4-11. Linearity Test of Nitrogen - Result Grid

Note. The linear regression of the $\delta^{15}\text{N}/^{14}\text{N}$ values vs. the working standard should be less than **0.06 ‰/V**.

Chapter 5

Carbon Measurement

5.1 Zero Enrichment of Carbon (Standard On/Off Test)

We assume that the user already has working experience with the ConFlo III Interface and IRMS. It is recommended to perform a simple check in order to test the analytical condition of ConFlo III and IRMS before measuring any samples. The most important checks to test the analytical condition are **zero enrichment** and **linearity test**.

1. Only as a guideline use the following Method from the File Browser's **Examples** folder (**CO2_zero.met**):

Instrument

The screenshot shows the 'Instrument' configuration window. The 'Reference Port' dropdown menu is set to 'Reference 1', which is highlighted by a red arrow. Other settings include 'Experiment' set to 'Continuous flow', 'Configuration' set to 'ConFlo', 'Gasconfiguration' set to 'CO2', and 'Acquisition Script' set to 'Acquisition.sct'. The 'Isotope MS' section contains 'Integration Time' at 0.200 [s], 'Peak Center Cup' at Cup 3, 'Peak Center Predelay (s)' at 20, and 'Peak Center Postdelay (s)' at 10. The 'Reference Device' section has 'Use Scripts' unchecked and 'Reference Port' set to Reference 1.

Figure 5-1. Zero Enrichment of Carbon - Instrument tab

- Select the reference port your reference gas is connected to at the ConFlo interface (e.g. Reference 1).
- Adapt the reference gas port to the respective column of the Time Events list: e.g. if you choose Reference 1, the on-off entries must occur in the Reference 1-On column of the Time Events list (see Figure 5-2).

Time Events

Time [s]	Start Sampler	Elemental Anal. - On	Dilution - On	Reference 1 - On	Reference 2 - On	Switch Gas
20				●		
40					●	
70				●		
90					●	
120				●		
140					●	
170				●		
190					●	
220				●		
240					●	
270				●		
290					●	
320				●		
340					●	
370				●		
390					●	

Figure 5-2. Zero Enrichment of Carbon - Time Events tab

- Recognize the eight on-off pulses shown above in the expected chromatogram (Figure 5-8).
- Note that the off-time of the e.g. fourth reference gas pulse is 190 s (Figure 5-4).

Acquisition Start	Immediately	Acquisition End Time [s]	420
-------------------	-------------	--------------------------	-----

Figure 5-3. Zero Enrichment of Carbon - Acquisition Start and Acquisition End Time

Evaluation

Instrument | Time Events | Evaluation@CO2 | Peak Detection@CO2 | Printout@CO2

Evaluation Type: >>

Ref. Nr.:	Ref. Time:	Ref. Name:	d 13C/12C	vs.	d 18O/16O	vs.
1	190.00	CO2_zero	0.000	VPDB	0.000	VSMOW

Reference/Blank

Significant Peak Start [s] Significant Peak Stop [s]

Amount Percent [%] Unit

Figure 5-4. Zero Enrichment of Carbon - Evaluation tab

Note. At **Ref. Time**, the **off**-value of an arbitrary reference gas peak according to the Time Events list must be typed (e.g. 190 s as the **off**-value of the forth reference gas peak, see Figure 5-2).

Peak Detection

Instrument | Time Events | Evaluation@CO2 | **Peak Detection@CO2** | Printout@CO2

Perform Peak Detection Perform Background Detection Detection on Mass

Detection Parameter

Start Slope [mV/s]

End Slope [mV/s]

Peak Min Height [mV]

Peak Resolution [%]

Max Peak Width [s]

Perform Timeshift

Background Parameter

Background Type

History [s]

Auto Square Pulse Recognition / Timeshift Suppression

Enable Factor rArea / Pk Width / Pk Height

Figure 5-5. Zero Enrichment of Carbon - Peak Detection tab

<< Advanced Parameter

Peak Detection Parameter Sets

Nr.:	Start Detection [s]	Stop Detection [s]
1	-1.000	-1.000

Smoothing

Smooth Type Number Of Datapoints

Figure 5-6. Zero Enrichment of Carbon - Advanced Parameters in Peak Detection tab

Note. A value of -1 denotes unlimited.

Printout

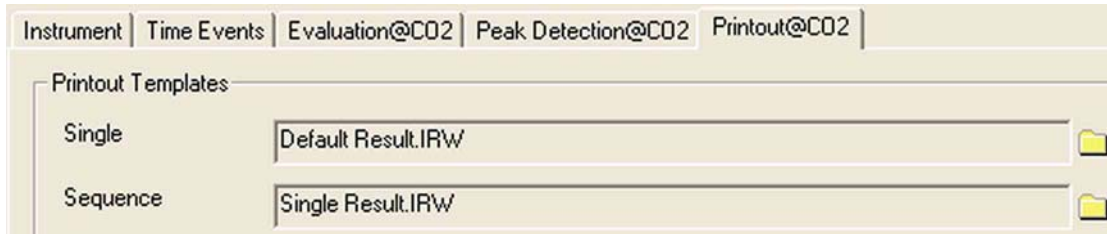


Figure 5-7. Zero Enrichment of Carbon - Printout tab

2. At the corresponding pressure regulator of ConFlo III, set the ion intensity of mass 44 (i.e. $^{12}\text{C}^{16}\text{O}^{16}\text{O}$) to 3 V - 4 V.
3. Create a new Sequence.
4. Press the **Start** button and confirm by **OK**.
5. Expected data after three or four measurements:

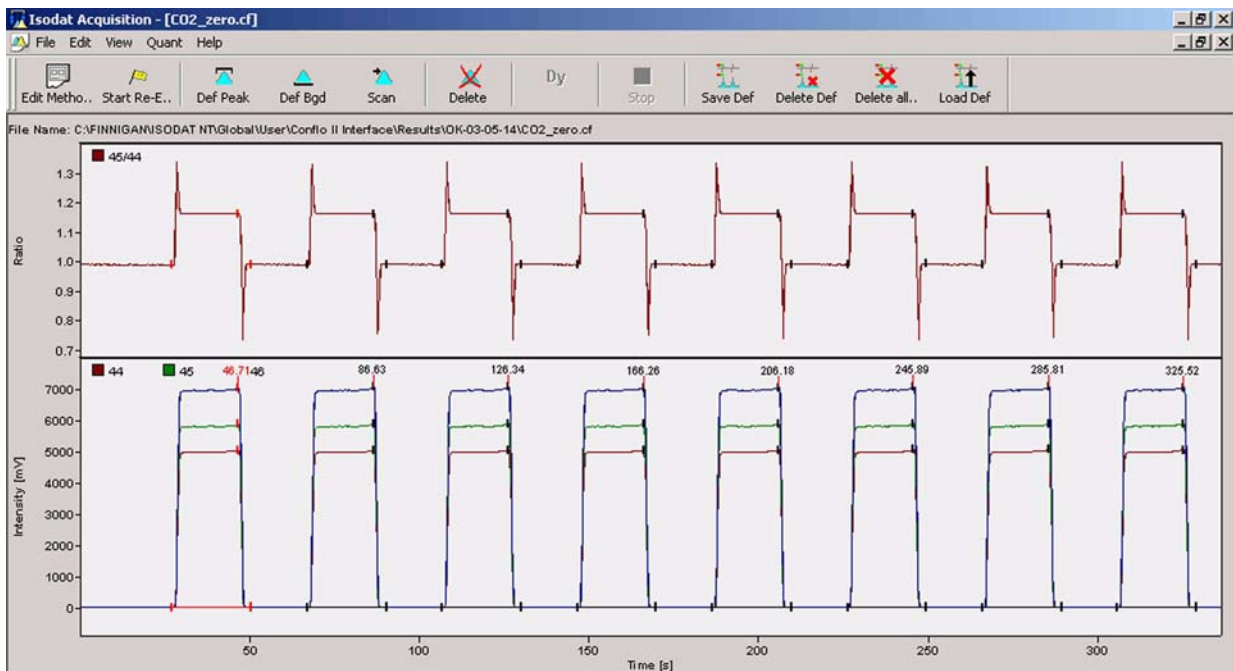


Figure 5-8. Zero Enrichment of Carbon - Chromatogram

Recognize the eight peaks shown above as the eight on-off pulses in the Time Events list (Figure 5-2).

CO2	Error	Extended	Sequence Line									
Peak Nr.	Start [s]	Rt [s]	Width [s]	Ampl. 44 [mV]	Ampl. 45 [mV]	BGD 44 [mV]	BGD 45 [mV]	Area All [Vs]	d 13C/12C [per mil] vs. VPDB	AT% 13C/12C [%]	d 18O/16O [per mil] vs. VSMOW	
1	27.3	46.7	23.2	5040	5875	3.9	2.7	96.047	0.055	1.105719	0.084	
2	67.0	86.6	23.4	5061	5897	4.0	2.8	96.074	0.006	1.105666	0.031	
3*	106.9	126.3	23.4	5061	5898	4.0	2.9	96.170	0.000	1.105659	0.000	
4	146.8	166.3	23.2	5053	5889	4.0	2.9	96.101	-0.019	1.105637	0.025	
5	186.5	206.2	23.4	5055	5890	4.0	2.9	96.144	-0.019	1.105638	-0.012	
6	226.5	245.9	23.2	5065	5904	4.0	2.9	96.209	-0.033	1.105622	-0.000	
7	266.2	285.8	23.4	5066	5903	4.0	2.9	96.132	-0.019	1.105638	0.004	
8	306.1	325.5	23.2	5056	5892	4.0	2.9	96.162	-0.006	1.105652	-0.001	

Figure 5-9. Zero Enrichment of Carbon - Result Grid

Note. The standard deviation of $\delta^{13}\text{C}/^{12}\text{C}$ should be **0.05 ‰** or less.


To obtain the standard deviation of all eight peaks perform as follows:

d 13C/12C [per mil] vs. VPDB
0.055
0.006
0.000
-0.019
-0.019
-0.033
-0.019
-0.006

- Click on the column header of the *d 13C/12C [per mil] vs. VPDB* column. It will be highlighted.



- Right-click on the column header.
- Choose *Calculate*.

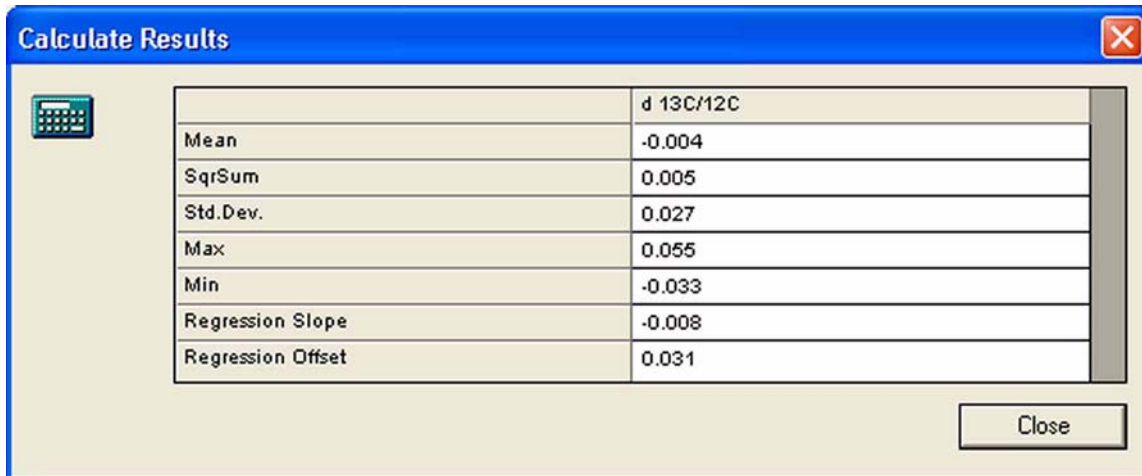

5.2 Linearity Test of Carbon

- Use the same method as defined for *zero enrichment* (i.e. *CO2_zero.met*, see chapter **Zero Enrichment of Carbon (Standard On/Off Test)** on page 5-2).
- Start the acquisition as single run.
- At each detection of a peak de- or increase the reference gas pressure at ConFlo III. Expected data after three or four measurements:

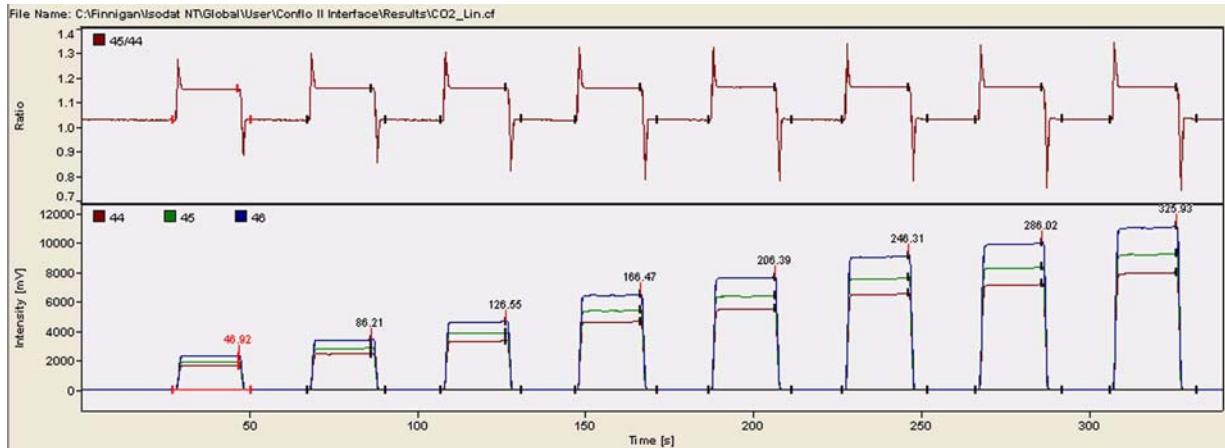


Figure 5-10. Linearity Test of Carbon - Chromatogram

CO2	Error	Extended	Sequence Line									
Peak Nr.	Start [s]	Rt [s]	Width [s]	Ampl. 44 [mV]	Ampl. 45 [mV]	BGD 44 [mV]	BGD 45 [mV]	Area All [Vs]	d 13C/12C [per mil] vs. VPDB	AT% 13C/12C [%]	d 18O/16O [per mil] vs. VSMOW	
1	27.5	46.9	23.0	1643	1910	12.3	15.4	31.347	-0.020	1.105637	0.126	
2	67.4	86.2	23.4	2450	2847	12.2	15.3	46.763	-0.002	1.105656	0.132	
3*	107.1	126.5	23.8	3343	3886	12.2	15.3	63.796	0.000	1.105659	0.000	
4	147.0	166.5	24.5	4660	5417	12.3	15.3	88.951	-0.035	1.105621	-0.123	
5	186.7	206.4	24.9	5514	6410	12.4	15.5	105.362	-0.083	1.105567	-0.213	
6	226.7	246.3	25.3	6583	7650	12.5	15.6	125.489	-0.123	1.105525	-0.349	
7	266.4	286.0	25.7	7203	8373	12.7	15.9	137.413	-0.181	1.105460	-0.422	
8	306.3	325.9	25.7	8011	9310	12.9	16.0	153.033	-0.223	1.105415	-0.580	

Figure 5-11. Linearity Test of Carbon - Result Grid

Note. The linear regression of the $\delta^{13}\text{C}/^{12}\text{C}$ values vs. the working standard should be less than **0.06 ‰/V**.

Chapter 6

Dual Measurement

6.1 Introduction

It is possible to perform dual measurements of nitrogen and carbon from a single sample with the system (ConFlo II/III and IRMS) in less than eight minutes offering fast sample throughput and high productivity. The technique is suitable for both solid and liquid samples.

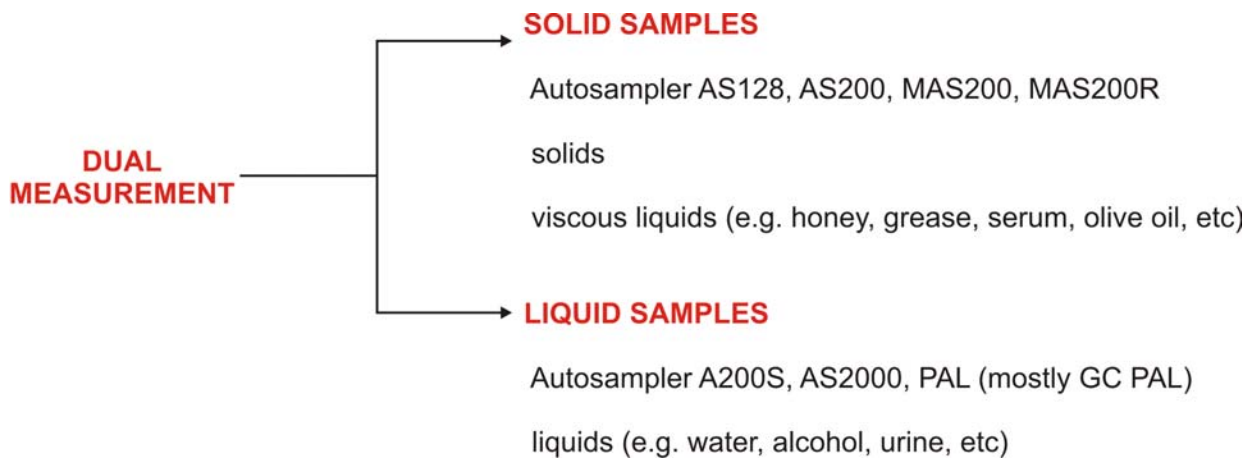


Figure 6-1. Schematic Diagram of Running Solid and Liquid Samples

As shown in Figure 6-1, different autosamplers are used to measure different sample types. Concerning the installation of the autosamplers refer to the manuals of the corresponding manufacturers.

6.2 Procedure

For the analysis of two isotopic species (nitrogen and carbon) from a single sample, a method, which comprises both gas configurations must be defined. The acquisition can be completed after less than eight minutes.

Due to the separation of the gases on a GC column, nitrogen appears first. As soon as the nitrogen peak (3) has been identified, Isodat 2.0 stops the nitrogen acquisition. The magnet “jumps” to the CO₂ configuration. A magnet jump means that the magnetic field is changed due to the entries from the mass calibration enabling the system to now collect masses 44, 45, 46 in the cups used for N₂ masses before.

If no nitrogen peak can be found, Isodat 2.0 waits a certain time, which has been set in the method, before automatically switching to CO₂ configuration.

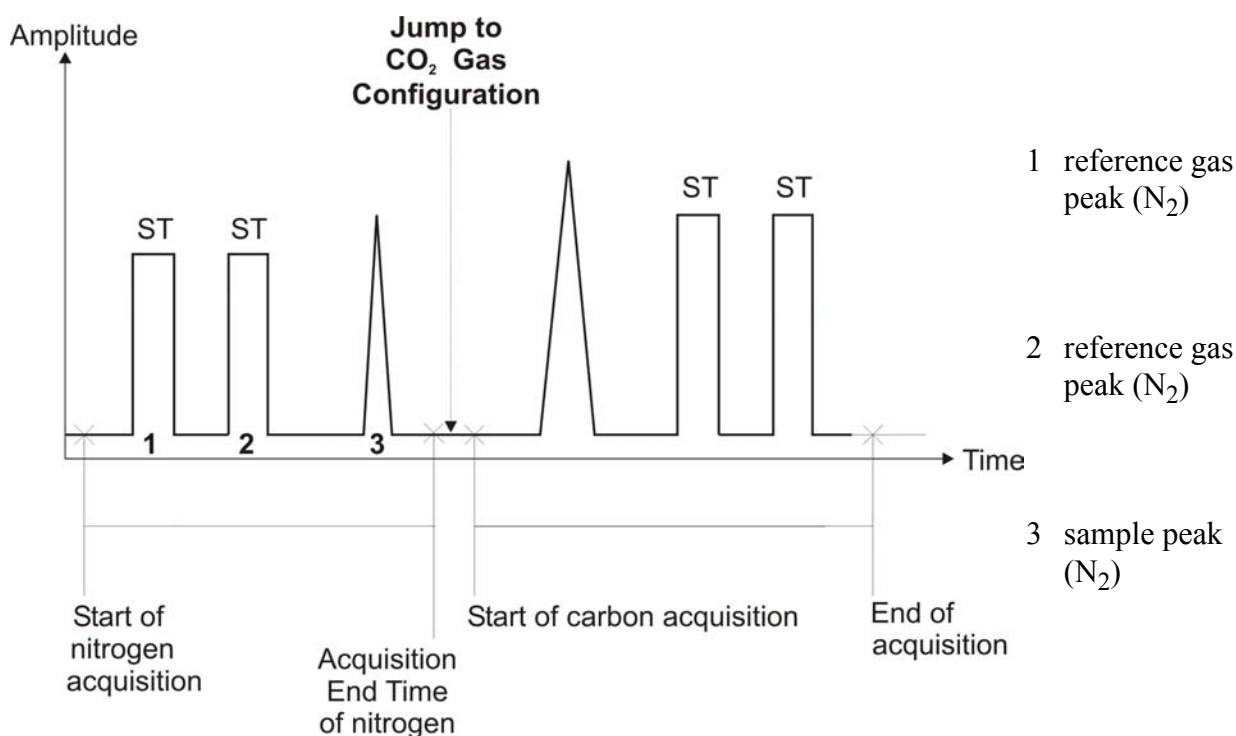


Figure 6-2. Dual Measurement - Schematic Chromatogram

6.3 How to Perform a Jump Calibration

To determine the isotope ratios of different elements during the same run, **switching to another Gas Configuration** is necessary.

In contrast to a single element measurement, in which the magnetic field runs the gamut from high to low and after that to the pre-calculated magnetic field, there is not sufficient time in Dual Measurement to perform this procedure for the next Gas Configuration.

For this reason, a **Jump Calibration** from the first Gas Configuration to the next Gas Configuration is necessary. It calibrates for a fast variation of the magnetic field.

After the Jump Calibration has been performed, the computer finds exactly the peak center even without performing any peak center procedure.

Temperature changes may influence the magnet performance on a daily basis. Therefore, a Jump Calibration should be performed daily.

Note. Both Gas Configurations used (i.e. N2 and CO2) need to be configured prior to Jump Calibration.

Note. A ConFlo configuration needs to be configured.



- Open **Instrument Control**.



- Switch to ConFlo **Configuration**.

The following window must appear at the **Accessories** toolbar. If it does not appear, check, if it is not visible (refer to Figure 6-3).

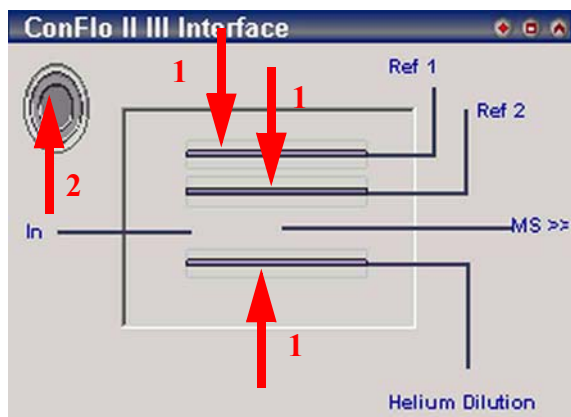


Figure 6-3. ConFlo II III Interface Window


The *ConFlo II III Interface* window allows:
 reference gas inlet and/or dilution (1)
 autosampler/Elemental Analyzer start signal (2)

Capillaries can be moved/activated by mouse clicks.

- Open **Reference 1** by a **click** on **Ref 1** capillary. If CO₂ is connected to **Reference 2**, click **Ref 2** capillary.

Thus, the blue capillary symbol elongates and CO₂ enters the system.

Comment: If the ConFlo II III Interface Window is not Visible

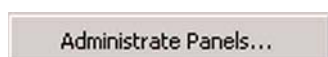

- **Right-click** somewhere into the *Accessories* toolbar.
- Click on the *Administrative Panels...* button.

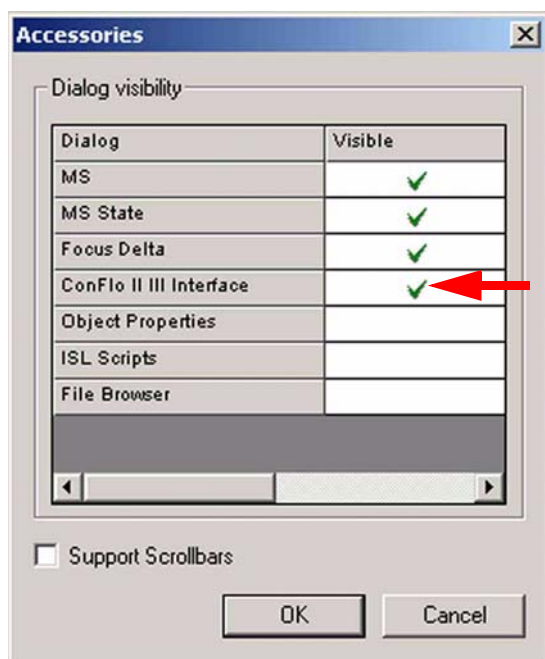


Figure 6-4. Visible Windows at the Accessories Toolbar


The *Accessories* window allows to mark dialogs (i.e. windows) by ✓. If you mark a dialog, the respective window will be shown at the *Accessories* toolbar.

- Mark *ConFlo II III Interface*.
- Confirm by **OK**.


Thus, the ConFlo II III Interface window will be shown at the *Accessories* toolbar.

Comment: If only the Title Bar of the ConFlo II III Interface Window is Visible



- Click on the *Unshrink* button  at the right corner of the Title bar.
Thus, ConFlo II III Interface window will appear.

Jump Calibration - Before Start

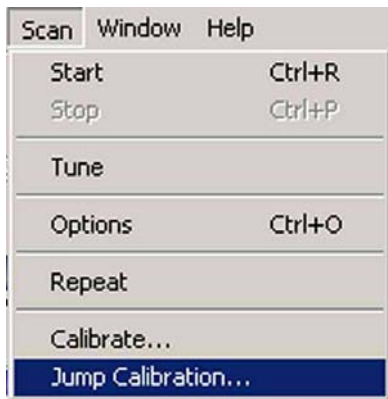

- Press *Scan > Jump Calibration...* or



- Alternatively, press the *Jump* button on the toolbar next to the Instrument Control window.

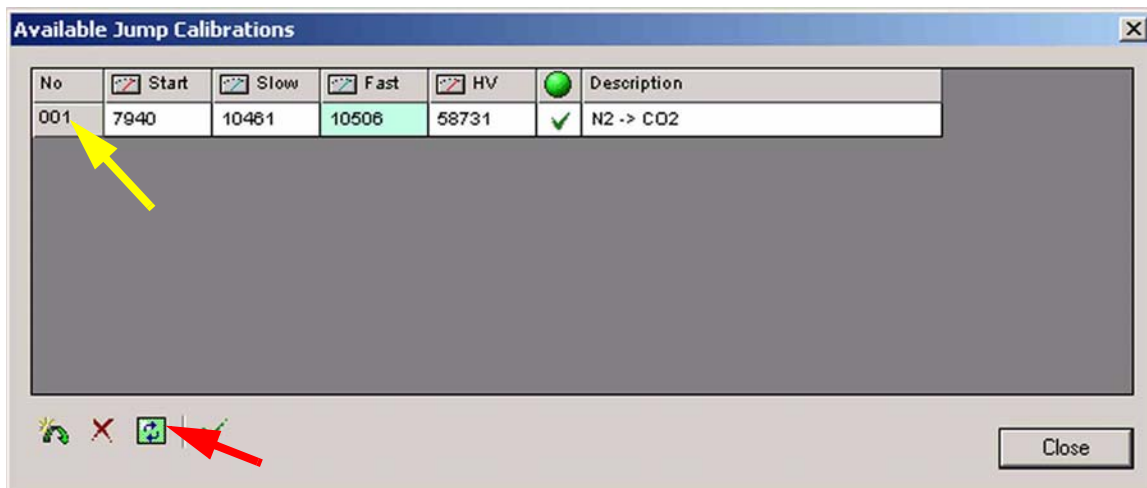


Figure 6-5. Available Jump Calibrations

- In the list of available Jump Calibrations mark the one for $N_2 \rightarrow CO_2$ by *clicking its No.*, e.g. 001.
- **Slow** compensate hysteresis by Max/Min settings of the magnet.
- **Fast** magnet setting of Jump Calibration
- **HV** high voltage setting of Jump Calibration

Note. Before pressing the **Recalibrate** button, make sure that CO_2 reference gas is activated, i.e. that CO_2 enters the source.



- Press the **Recalibrate** button.

Comment: If no Jump Calibration is Available

If *no* Jump Calibration is available in the list, *create a new one* as follows:

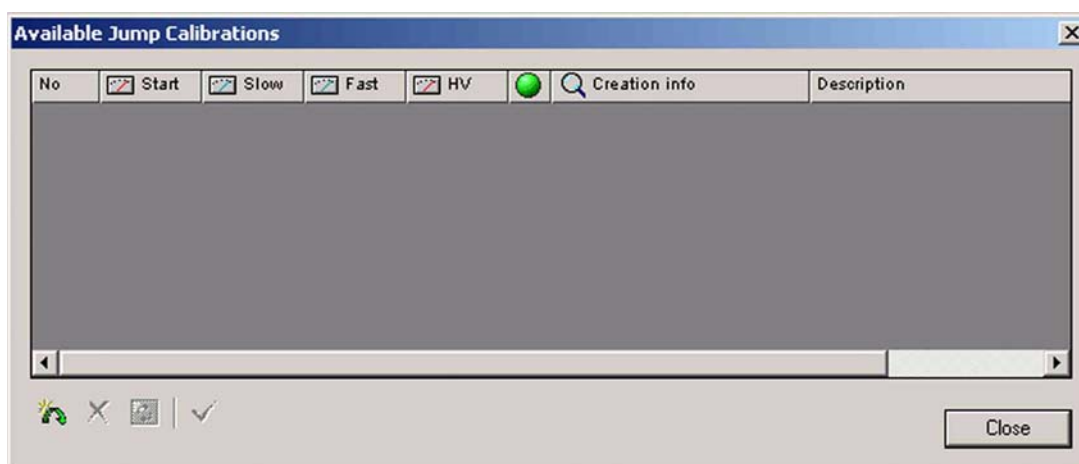


Figure 6-6. No Jump Calibration Available



- Press the **New** button.



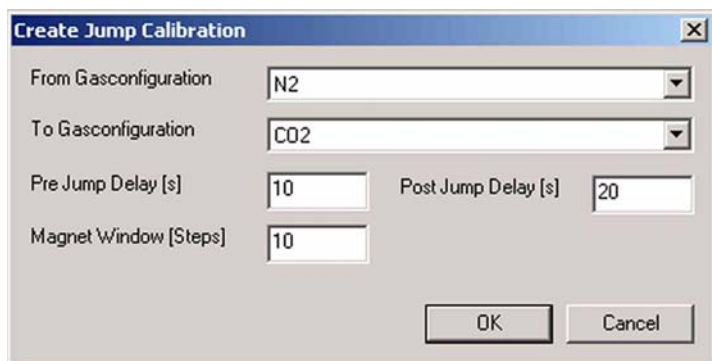


Figure 6-7. Define the Parameters of the new Jump Calibration

- At *From Gas Configuration* select *N2*.
- At *To Gas Configuration* select *CO2*.
- At *Pre Jump Delay [s]* accept the default value.
- At *Post Jump Delay [s]* accept the default value.
- At *Magnet Window [Steps]* accept the default value.
- Confirm by *OK*.

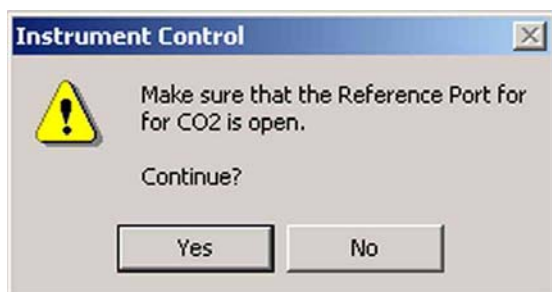



- Confirm by *Yes*, if you have already opened the Reference Port for CO₂ as shown in Figure 6-3.

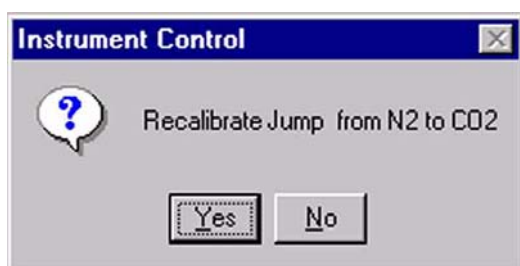

- Confirm by *Yes* to start Jump Calibration.

Jump Calibration - Procedure

- 1a Jump to **CO₂** (along hysteresis curve).
- 1b Perform a peak center for **CO₂** to get the signal height.
- 2 Jump to **N₂** (along hysteresis curve; **N₂** is origin).
- 3a Jump to **CO₂** (*not* along hysteresis curve).
- 3b Perform a peak center for **CO₂** to catch the peak.

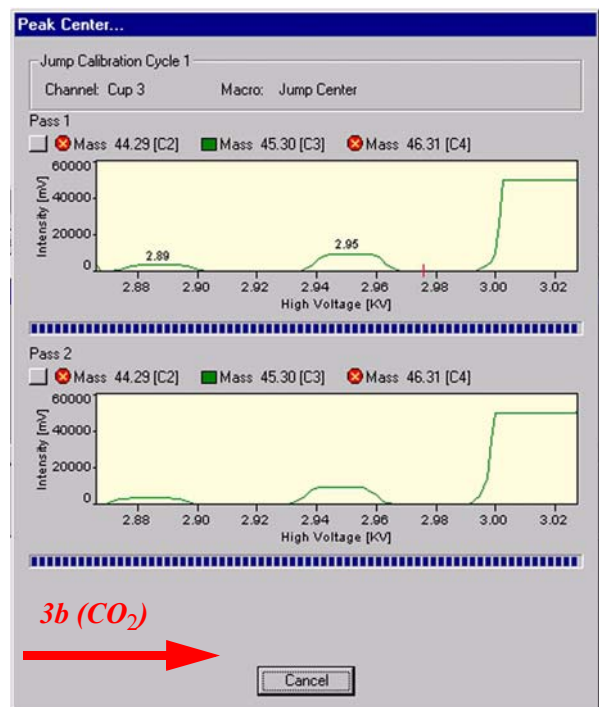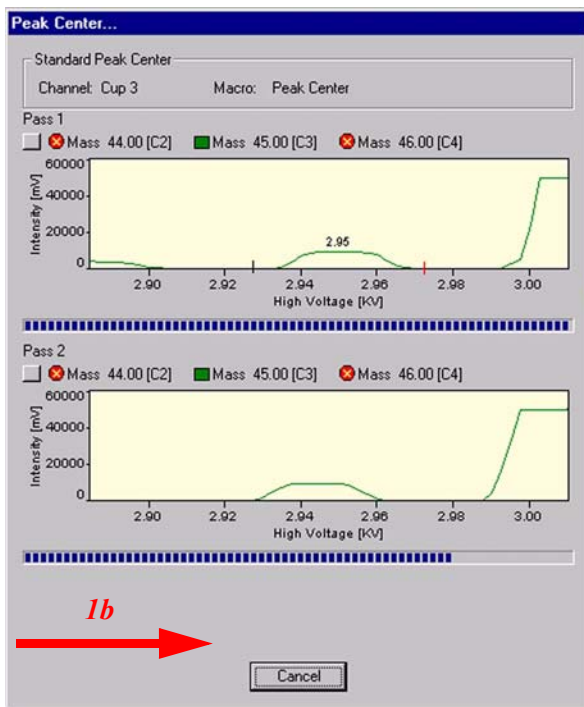

Chapter 7

Sulfur Measurement

7.1 Introduction

In comparison to isotope ratio measurements of nitrogen (^{15}N) and carbon (^{13}C) in organic and inorganic matter, the analysis of sulfur (^{34}S) has always been difficult. Analyzing biological sulfur makes experimental difficulties caused by the low abundance of sulfur in organisms (e.g. 0.2 wt % [mg/mg] in plants), as well as by the fact that sulfur is present as a mixture of organic and inorganic compounds. Difficulties in using the Elemental Analyzer arise due to a large amount of carbon in the same samples, which quickly exceeds the capacity of combustion reactors. This causes incomplete combustion.

Thermo Electron (Bremen) has developed a technique for precise, accurate and fast sulfur measurement, which puts it on a par with carbon and nitrogen in terms of ease of use and sample size.




Due to the high natural abundance of the heavier isotope ^{34}S , less amplification is required and may become necessary for all Delta IRMS before Delta^{plus} XP (since 2002). It is recommended to use a smaller resistor ($1 * 10^{10} \Omega$) on the cup for mass 66 (usually $3 * 10^{10} \Omega$).

7.2 Procedure

Sulfur measurements are performed using a specially equipped Elemental Analyzer. Combustion and reduction are carried out in a single reactor filled with tungsten oxide (WO_3) and copper (Cu) as reducing agent. Alternatively, CuO can be used instead of WO_3 .

The technique used for sulfur determination is based on the quantitative "Dynamic Flash Combustion" method. The samples - sometimes together with V_2O_5 - are wrapped in tin capsules and placed into the autosampler. Then they are continuously purged with helium (or oxygen using NA 2500 or EA 1110) to remove any traces of water and nitrogen. When a sample is dropped into the reactor, the helium stream is temporarily enriched with pure oxygen. The sample and its container melt as the tin promotes a violent reaction (flash combustion). Under these favorable conditions, even thermally resistant substances are completely oxidized.

In the reactor, e.g. BaSO_4 is thermally decomposed within a tin capsule. The following reactions can then take place (see Bailey, S.A. and Smith, J.W., 1972):

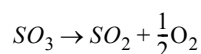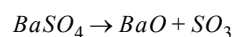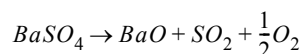

Note. Bailey, S.A. and Smith, J.W. (1972): An improved method for the preparation of sulfur dioxide from barium sulfate for isotope ratio studies. *Anal. Chem.* **44**, 1542-1543.

Although the process does not require oxygen, better combustion has been experienced when O_2 is injected and vanadium pentoxide (V_2O_5) is added to the sample. If either the O_2 pressure is low or a bad catalyst is selected or the reactor has too much ash, combustion will proceed slowly. A slow stream of SO_2 through the system causes adsorption at the tubing wall.

Note. It is important that enough reduced copper is present in the combustion tube. If this is not assured, SO_3 will only be reduced partially and isotope fractionation will occur.

"Light" isotopes ($^{32}\text{SO}_3$) are reduced more easily than "heavy" isotopes ($^{34}\text{SO}_3$). Therefore, SO_2 gas is depleted in ^{34}S compared to the original sample, and the δ value becomes more negative.

7.3 Sulfur Measurement Kits

To perform a sulfur measurement you can get two *sulfur measurement kits*, provided by Thermo Electron (Bremen):

- Sulfur Measurement Kit for ConFlo III
- Sulfur Measurement Kit for Flash Elemental Analyzer

Sulfur Measurement Kit for ConFlo III

It consists of the following parts:

Table 7-1. Parts of Sulfur Measurement Kit for ConFlo III (Part No. 115 7100)

Quantity	Description	Part No.
1	attachment for exhaust tube	112 1390
1	self-adhesive heating foil for ConFlo III	114 1180
1	power supply for self-adhesive heating foil for ConFlo III	204 8580

Sulfur Measurement Kit for Flash Elemental Analyzer

It consists of the following parts:

Table 7-2. Parts of Sulfur Measurement Kit for Flash Elemental Analyzer (Part No. 115 7110)

Quantity	Description	Part No.
1	sulfanilamide	106 9140
1	Teflon GC column	114 1170
5	Al olive (2 mm)	114 1210
5	nut	112 1370
3.5 m	Teflon tube	114 1220*
2	combustion reactor (packed, "ready for use")	111 8121
3	union	114 1230
0.05 m	capillary tube (1/16" * 0.8 mm)	060 5470

*For additional part numbers (e.g. of nuts) see **Installing the Teflon Tubing** on page 7-6 and especially Figure 7-2.



Warning. When working with sulfur dioxide (SO₂), good ventilation is essential. Otherwise, the gas can be hazardous to your health.



Warning. To ensure operational safety, a CO, SO₂ and H₂ detector with an alarm must be installed!

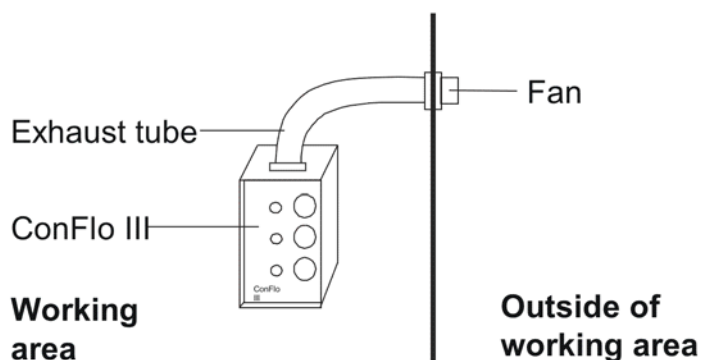


Figure 7-1. ConFlo III Interface with Exhaust Tube and Fan

The exhaust tube should be installed on top of the ConFlo III interface as shown in Figure 7-1 to remove the toxic sulfur dioxide (SO₂) from inside of ConFlo III out of your working area.

7.4 Preparing the System for a Sulfur Measurement

1. Install the exhaust tube on ConFlo III as shown in Figure 7-1.
2. Install the self-adhesive heating foil as shown in Figure 7-3.
3. Install the properly packed reactor within the Elemental Analyzer.
4. Replace all stainless steel tubing with Teflon tubing as shown in Figure 7-2.
5. Install the SO₂ GC column.

Installing the Teflon Tubing

To perform a sulfur measurement, a part of the stainless steel tubing needs to be replaced by Teflon tubing. The scheme below provides an overview:

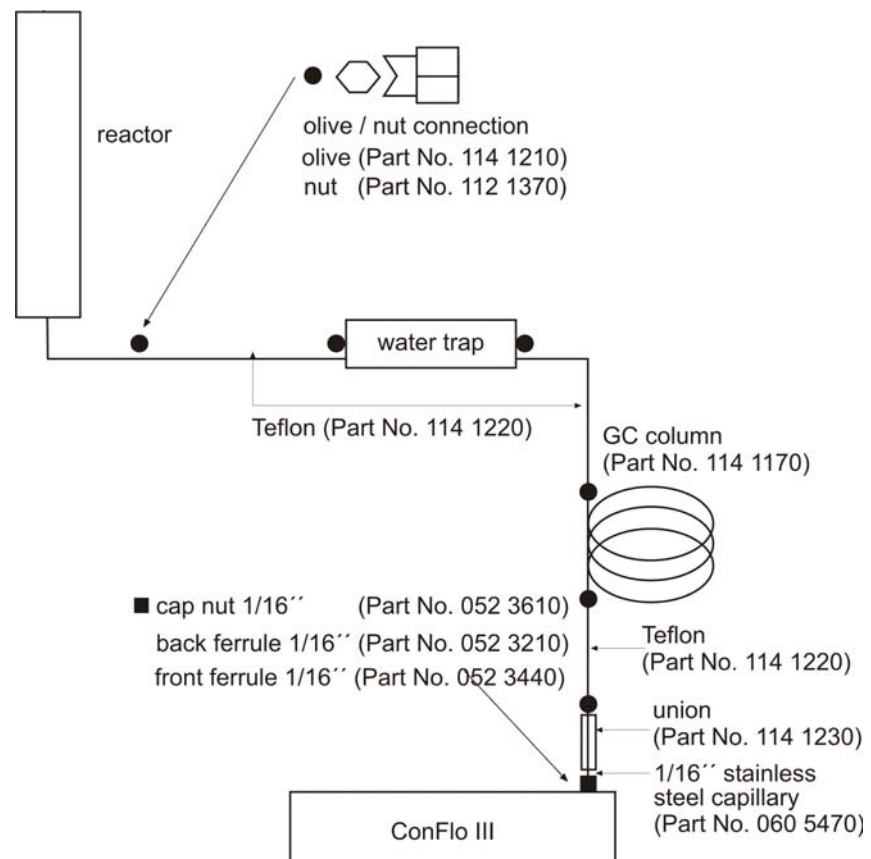


Figure 7-2. Installing the Teflon Tubing

Installing the Self-Adhesive Heating Foil

Note. SO₂ is liquid at higher pressures. Therefore, the manometer should be heated to avoid condensation of SO₂. The temperature must be between 60 °C and 70 °C. It may vary with ventilation.

1. Remove the backing paper from the self-adhesive heating foil.
2. Paste the heating foil in the middle of the manometer's rear side (of Ref 2) as shown in Figure 7-3.

Note. Take care that the wires of the heating foil point downwards when pasting it on the manometer's rear side.

3. Simply insert the wires into the plug socket where they fix themselves.


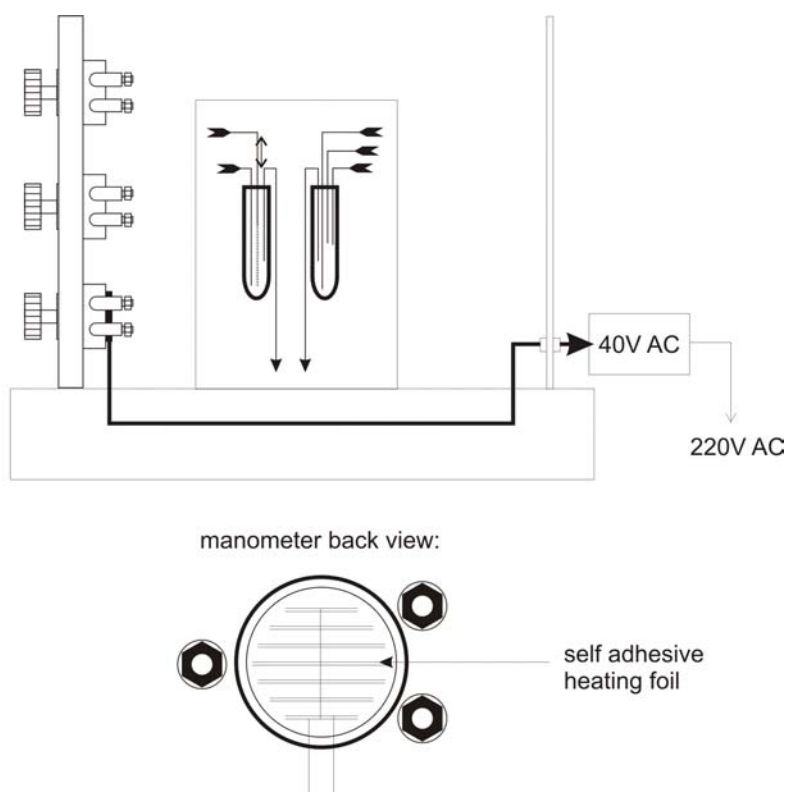


Figure 7-3. Installing the Self-Adhesive Heating Foil

Reactor Filling

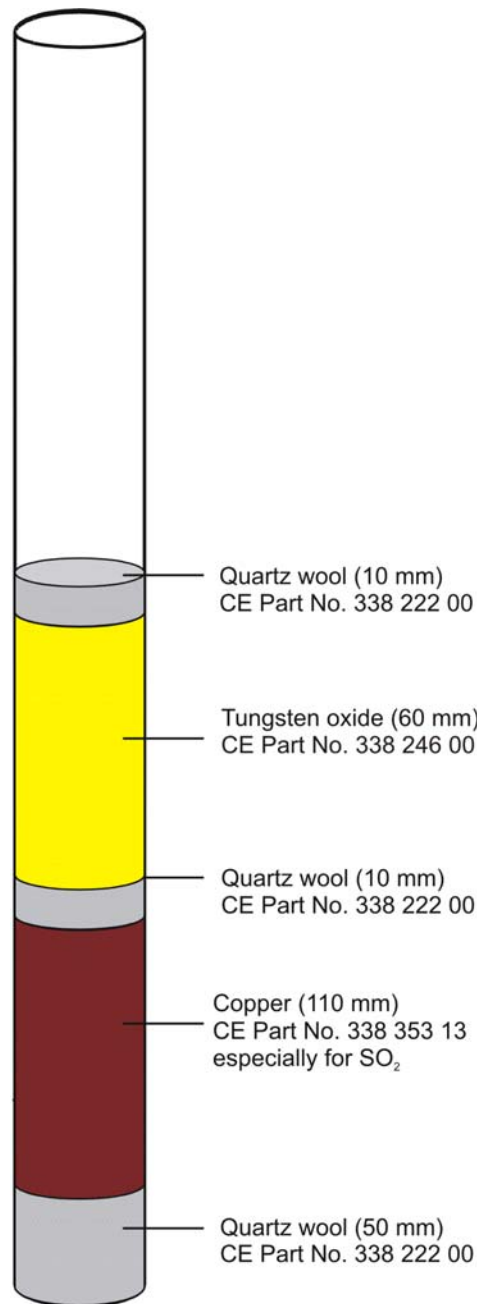

We recommend to use a Thermo Electron "ready for use" reactor (Part No. 111 8121). The reactor is 470 mm long. If the reactor must be packed by the user, the packing should be as shown left.

About 150 to 200 analyses can be performed using the reactor type described above. With the exception of tungsten oxide, no parts of the reactor can be used a second time.

Figure 7-4. Packing of the Reactor

***Option.** Instead of tungsten oxide also 60 mm of copper oxide (CE Part No. 338 217 10) can be used for the packing of the reactor.

7.5 Before Starting a Sulfur Measurement

We assume that the user already has working experience with the isotope ratio mass spectrometer. Before starting a sulfur measurement make sure that:

1. The Elemental Analyzer is set up, that is:
 - The SO₂ reactor is packed and installed.
 - The water trap is installed.
2. The special SO₂ stainless steel (or Teflon) GC column is installed.
3. The connection between reactor, water trap, GC column and ConFlo III is made up of Teflon tubing.
4. The gases needed (He, O₂) are available and have been connected to the corresponding positions.
5. Leak-check all connections outside of ConFlo III by brushing all fittings carefully with soapsuds.
6. The exhaust tube is connected to the ConFlo III and ventilation is working.
7. SO₂ is connected to Ref 2 of ConFlo III.
8. The self-adhesive heating foil is installed / switched on.
9. Compressed air is connected to ConFlo III and Elemental Analyzer.
10. ConFlo III configuration is installed and has already been in use for e.g. nitrogen or carbon analysis.
11. The IRMS has been calibrated for SO₂ measurements.
12. Switch the Elemental Analyzer on.
13. Adjust the gas pressure of Elemental Analyzer and ConFlo III
Take the following settings as a guideline.

For EA 1108, EA 1110 and NA 2500

Table 7-3. Settings for EA 1108, EA 1110 and EA 2500

He	Elemental Analyzer	100 kPa
O ₂	Elemental Analyzer	150 kPa
Purge with	O ₂	40 ml/min
Purge with	He	110 ml/min (depending on the autosampler)
He	ConFlo III	1 bar
Ref. 2	ConFlo III	1 bar

For Flash Elemental Analyzer

Table 7-4. Settings for Flash Elemental Analyzer in Eager Software

Carrier	300 ml/min
Oxygen	175 ml/min
Reference	110 ml/min (depending on autosampler)
Cycle run time	60 s
Sampling delay	23 s
Oxygen Injection End	3 s

14. Switch the ion source **OFF** and inject the He-dilution capillary.
15. Slowly open the needle valve. Make sure that the needle valve is heated.
16. Make sure that the ion source heater and the change over valve heater are operating, if available.
17. Switch the ion source **ON**.
18. After approximately 30 min. the expected background at the respective reactor and column temperature should be as follows:

Stainless Steel GC Column

Table 7-5. Expected Background Using a Stainless Steel GC Column

reactor temperature	column temperature	mass cup range	28 mV	29 mV	Ar mV	H ₂ O V	64 mV	66 mV
30 °C	30 °C		50	60	180	1.5		
400 °C	100 °C		110	140	350	1.8		
1020 °C	100 °C		70	65	220	1.2	< 5	< 10

Note. Background values may vary depending on sensitivity and focus settings. They are given here as a general guideline.

Teflon GC Column

Table 7-6. Expected Background Using a Teflon GC Column

reactor	column	mass	28	29	Ar	H ₂ O	64	66
temperature	temperature	cup	28	29	29	29	64	*66
		range	mV	mV	V	V	mV	mV
30 °C	30 °C		295	223	0.93	3.9		
400 °C	100 °C		422	337	1.30	5.6		
1020° C	100 °C		330	243	0.99	2.1	<5	<10

(* R_i=10¹⁰ Ω)

Note. Background values may vary depending on sensitivity and focus settings. They are given here as a general guideline.

Note. Teflon tubes are not absolutely tight against atmosphere. Therefore, the backgrounds of argon and nitrogen are higher than those of carbon and nitrogen measurements. Nevertheless, take care of leaks.

Note. When using normal amplification (i.e. resistor of 3 * 10¹⁰ Ω on cup for mass 66), the background values will be three times higher.

7.6 Create a Gas Configuration for a Sulfur Measurement

A Gas Configuration determines a combination of masses, which are collected in the cups, for evaluation of ratios and eventually δ values. The Gas Configuration is specific for the particular gas and is combined with a magnet field value taken from the mass calibration of your IRMS. The ratio groups determine the reported ratios of predefined masses.

Prior to defining this Gas Configuration ensure that the connected IRMS has the cups for the simultaneous detection of masses 64 and 66 and mass calibration for these cups has already been performed.


For a ^{34}S measurement, a Gas Configuration must be available for the masses 64 (i.e. $^{34}\text{S}^{16}\text{O}^{16}\text{O}$) and 66 (i.e. $^{32}\text{S}^{16}\text{O}^{16}\text{O}$). Otherwise, it must be created as follows.



- Open the *Acquisition* module.



- Open the *Gas Configuration Editor*. It is only available if no acquisition is running.

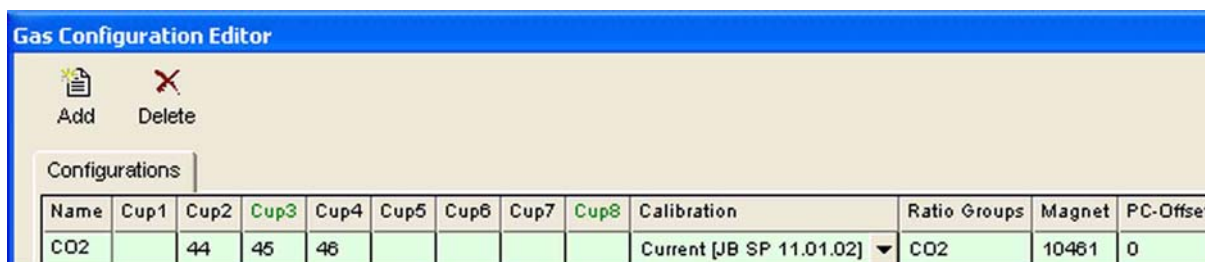

- Per default, the Gas Configuration *CO2* is being created as the first one.
- If the Gas Configuration *SO2* has already been created, it occurs in the list above.
- If the Gas Configuration *SO2* has not been created yet, it does not occur in the list above. Then, proceed as follows.



New Gasconfiguration

Name:

Template:

- **Add** a new Gas Configuration.
- Type **SO2** for the **Name**.
- Select a Gas Configuration as **Template** (e.g. **CO2**).

In the pulldown menu, only the already existing Gas Configurations are displayed. When creating the first Gas Configuration, **CO2** is displayed.

- Confirm by **OK**.

Isodat Acquisition

Confirm Formula.

Name : <SO2>
Formula: <CO2>

Press <Yes> for continue or <No> for edit Formula.

- Type **No**.

If you would type **Yes**, this would automatically mark the template (i.e. **CO2**) instead of **SO2** in the **Ratio Groups** window below.

Ratio Groups

Evaluations	Enable
CO2	<input type="checkbox"/>
CO	<input type="checkbox"/>
H2	<input type="checkbox"/>
N2	<input type="checkbox"/>
O2	<input type="checkbox"/>
N2O	<input type="checkbox"/>
SO2	<input checked="" type="checkbox"/>
CH3Cl	<input type="checkbox"/>
SF6	<input type="checkbox"/>
Argon	<input type="checkbox"/>
Air	<input type="checkbox"/>
SO-SO2	<input type="checkbox"/>
SE5	<input type="checkbox"/>

- Mark **SO2**.
- If Ratio Groups other than **SO2** are marked, unmark them all.

- Confirm by **OK**.


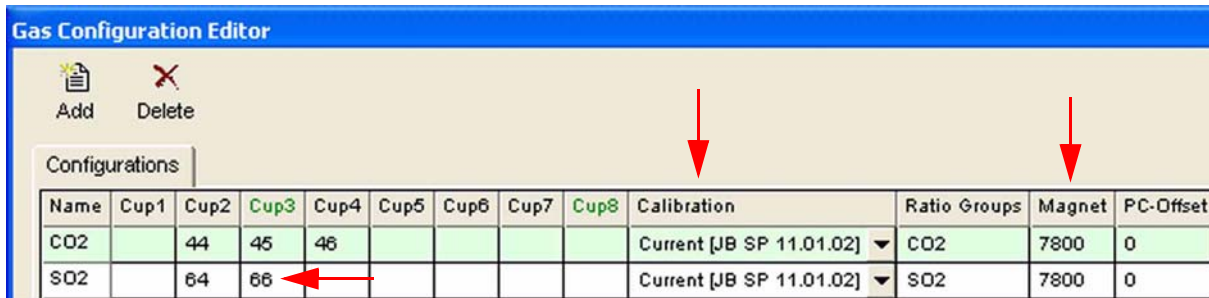


Figure 7-5. Creating a New Gas Configuration


- The new Gas Configuration **SO2** appears in the list as a row of its own.
- In the **Calibration** column select your current calibration file.

Note. The above figure shows a common cup configuration as used in most Delta mass spectrometers, i.e. universal triple collector.
If you have a special cup configuration, the respective masses will be collected in other cups!


- Fill in the correct masses (64 and 66 replace e.g. 44, 45 and 46) to the appropriate cups specific for your IRMS.

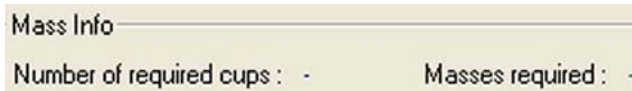

When highlighting the specific gas configuration by a click on its row, the number of cups required for measurement is displayed together with the assigned masses.

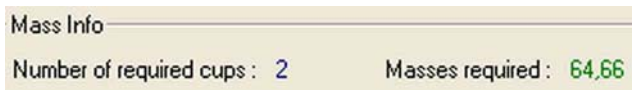

- Select a **Calibration**, which is valid for the selected cups.

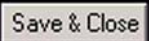Press the **Save & Close** button  .

7.7 Zero Enrichment of Sulfur (Standard On/Off Test)

We assume that the user already has working experience with the ConFlo III Interface and IRMS. It is recommended to perform a simple check in order to test the analytical condition of ConFlo III and IRMS before measuring any samples. The most important checks to test the analytical condition are *zero enrichment* and *linearity test*.

1. Only as a guideline use the following method from the File Browser's *Examples* folder (*SO2_zero.met*):

Instrument

The screenshot shows the 'Instrument' tab of the software interface. The 'Reference Port' dropdown menu is set to 'Reference 2', which is highlighted by a red arrow. Other visible settings include 'Experiment' set to 'Continuous flow', 'Configuration' set to 'ConFlo', 'Acquisition Script' set to 'Acquisition.sct', and 'Isotope MS' settings for 'Integration Time' (0.200 [s]), 'Peak Center Cup' (Cup 3), 'Peak Center Predelay (s)' (20), and 'Peak Center Postdelay (s)' (10).

Figure 7-6. Zero Enrichment of Sulfur - Instrument Tab

- Select the reference port your reference gas is connected to at the ConFlo interface (e.g. Reference 2).
- Adapt the reference gas port to the respective column of the Time Events list: e.g. if you choose Reference 2, the on-off entries must occur in the Reference 2-On column of the Time Events list (see Figure 7-7).

Time Events

Time [s]	Start Sampler	Elemental Anal. - On	Dilution - On	Reference 1 - On	Reference 2 - On	Switch Gas
40					●	
60						●
110					●	
130						●
180					●	
200						●
250					●	
270						●
320					●	
340						●
390					●	
410						●
460					●	
480						●
530					●	
550						●
600					●	
620						●

Figure 7-7. Zero Enrichment of Sulfur - Time Events Tab

Acquisition Start Acquisition End Time [s]

Figure 7-8. Zero Enrichment of Sulfur - Acquisition Start and Acquisition End Time

- Recognize the nine on-off pulses shown above in the expected chromatogram (Figure 7-12).
- Note that the off-time of the e.g. fourth reference gas pulse is 270 s (Figure 7-9).

Evaluation

Instrument | Time Events | Evaluation@SO2 | Peak Detection@SO2 | Printout@SO2

Evaluation Type: >>

Ref. Nr.:	Ref. Time:	Ref. Name:	d 18O/16O	vs.	d 34S/32S	vs.
1	270.00	SO2_zero	0.000	VSMQW	0.000	VCDT

Reference/Blank

Significant Peak Start [s] Significant Peak Stop [s]

Amount Percent [%] Unit

Figure 7-9. Zero Enrichment of Sulfur - Evaluation Tab

Note. At **Ref. Time**, the **off**-value of an arbitrary reference gas peak according to the Time Events list must be typed (e.g. 270 s as the **off**-value of the forth reference gas peak, see Figure 7-7).

Peak Detection

Instrument | Time Events | Evaluation@SO2 | Peak Detection@SO2 | Printout@SO2

Perform Peak Detection Perform Background Detection Detection on Mass

Detection Parameter

Start Slope [mV/s]

End Slope [mV/s]

Peak Min Height [mV]

Peak Resolution [%]

Max Peak Width [s]

Perform Timeshift

Background Parameter

Background Type

History [s]

Auto Square Pulse Recognition / Timeshift Suppression

Enable Factor rArea / Pk Width / Pk Height

<< Advanced Parameter

Peak Detection Parameter Sets

Nr.:	Start Detection [s]	Stop Detection [s]
1	-1.000	-1.000

Smoothing

Smooth Type

Number Of Datapoints

Figure 7-10. Zero Enrichment of Sulfur - Peak Detection Tab

Note. A value of -1 denotes unlimited.

Printout

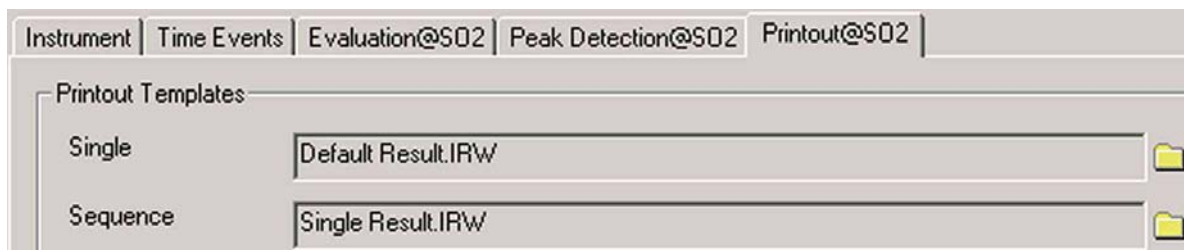


Figure 7-11. Zero Enrichment of Sulfur - Printout Tab

2. At the corresponding pressure regulator of ConFlo III, set the ion intensity of mass 64 (i.e. $^{32}\text{S}^{16}\text{O}^{16}\text{O}$) to approximately 3 V.
3. Create a new Sequence.
4. Press the **Start** button and confirm by **OK**.
5. Expected data after three or four measurements:

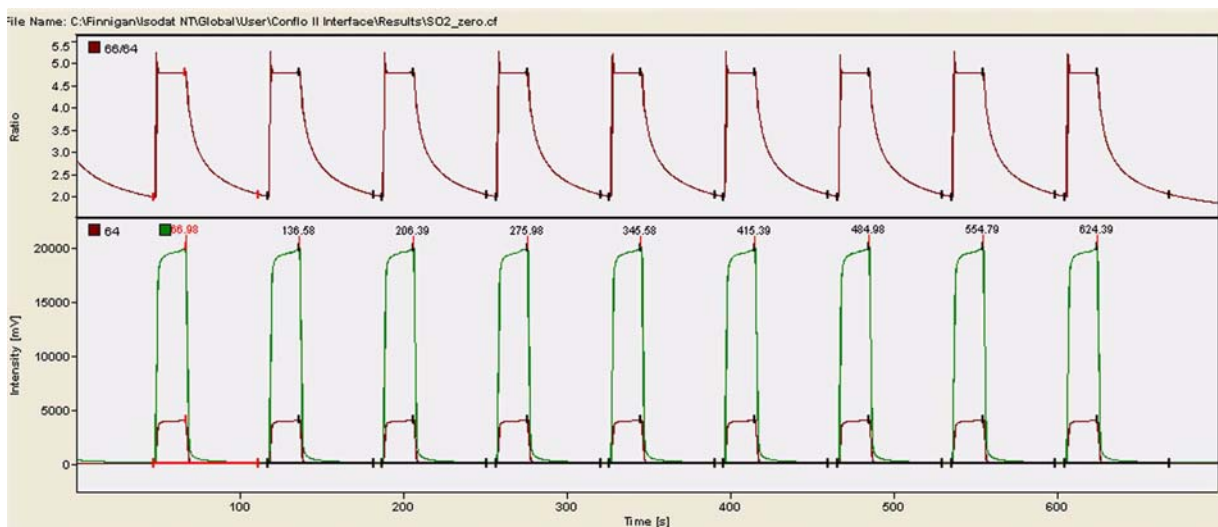


Figure 7-12. Zero Enrichment of Sulfur - Chromatogram

Recognize the nine peaks shown above as the nine on-off pulses in the Time Events list (Figure 7-7).

SO2		Error	Extended	Sequence Line								
Peak Nr.	Start [s]	Rt [s]	Width [s]	Ampl. 64 [mV]	Ampl. 66 [mV]	BGD 64 [mV]	BGD 66 [mV]	Area All [Vs]	Amt% [%]	d 66SO2/64SO2 [per mil] vs. SO2_zero	d 34S/32S [per mil] vs. VCDT	
1	47.8	67.0	63.7	4085	19953	33.5	163.9	79.345	-	0.021	0.022	
2*	117.4	136.6	64.4	4061	19838	34.1	166.8	79.141	-	0.000	0.000	
3	187.2	206.4	64.0	4070	19878	34.0	166.3	79.308	-	-0.040	-0.044	
4	256.8	276.0	64.4	4072	19887	33.9	165.8	79.329	-	-0.042	-0.046	
5	326.4	345.6	64.2	4075	19904	33.9	165.9	79.465	-	-0.048	-0.052	
6	396.2	415.4	63.7	4068	19869	33.8	165.8	79.164	-	-0.076	-0.083	
7	465.8	485.0	63.7	4094	19995	33.8	165.8	79.399	-	-0.078	-0.085	
8	535.6	554.8	63.7	4091	19981	33.8	165.9	79.386	-	-0.105	-0.114	
9	605.2	624.4	64.0	4083	19944	33.9	166.2	79.401	-	-0.117	-0.128	

Figure 7-13. Zero Enrichment of Sulfur - Result Grid

Note. The standard deviation of $\delta^{66}\text{SO}_2/^{64}\text{SO}_2$ should be 0.05 ‰ or less.

To obtain the standard deviation of all nine peaks perform as follows:

d 66SO2/64SO2
0.021
0.000
-0.040
-0.042
-0.048
-0.076
-0.078
-0.105
-0.117

- Click on the column header of the *d 66SO2/64SO2 [per mil] vs. SO2_zero* column. It will be highlighted.



- Right-click on the column header.
- Choose *Calculate*.

A screenshot of the 'Calculate Results' dialog box. It contains a table with statistical data for the column 'd 66SO2/64SO2'. A 'Close' button is located at the bottom right.

	d 66SO2/64SO2
Mean	-0.054
SqrSum	0.017
Std.Dev.	0.045
Max	0.021
Min	-0.117
Regression Slope	-0.016
Regression Offset	0.028

7.8 Linearity Test of Sulfur

- Use the same method as defined for *zero enrichment*, e.g. *SO2_zero.met*. See **Zero Enrichment of Sulfur (Standard On/Off Test)** on page 7-15.
- Start the acquisition as a single run.
- At each detection of a peak de- or increase the reference gas pressure at ConFlo III.

Expected data after three or four measurements:

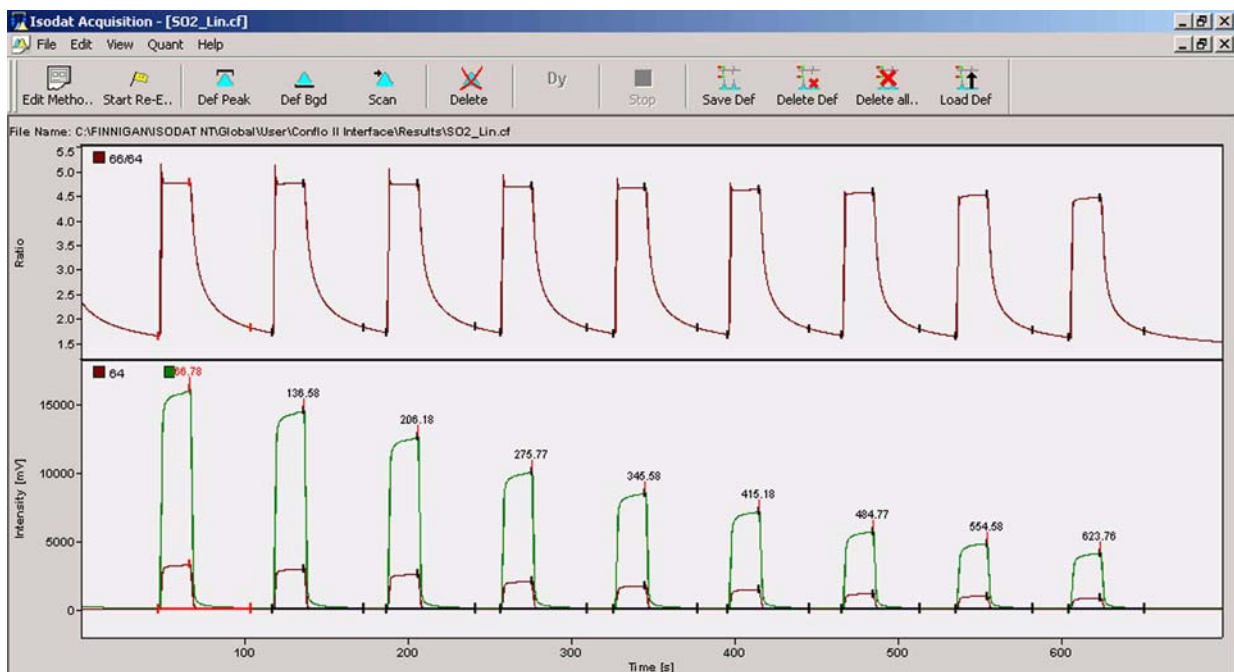


Figure 7-14. Linearity Test of Sulfur - Chromatogram

SO2	Error	Extended	Sequence Line								
Peak Nr.	Start [s]	Rt [s]	Width [s]	Ampl. 64 [mV]	Ampl. 66 [mV]	BGD 64 [mV]	BGD 66 [mV]	Area All [Vs]	Amt% [%]	d 66SO2/64SO2 [per mil] vs. SO2_zero	d 34S/32S [per mil] vs. VCDT
1	47.5	66.8	56.6	3289	16053	22.7	102.8	63.885	-	-0.008	-0.009
2	117.4	136.6	56.0	2964	14464	24.8	113.7	57.796	-	-0.031	-0.034
3*	186.9	206.2	54.5	2579	12586	25.2	115.4	50.025	-	0.000	0.000
4	256.8	275.8	53.1	2046	9986	24.9	114.2	39.707	-	0.038	0.042
5	326.4	345.6	51.2	1728	8433	24.2	110.8	33.495	-	0.099	0.108
6	396.2	415.2	49.5	1448	7066	23.6	107.6	28.036	-	0.131	0.142
7	465.8	484.8	48.1	1145	5590	22.9	104.1	22.150	-	0.260	0.283
8	535.6	554.6	47.2	972	4744	22.1	100.0	18.758	-	0.288	0.314
9	605.2	623.8	45.8	827	4034	21.5	96.8	15.917	-	0.359	0.392

Figure 7-15. Linearity Test of Sulfur - Result Grid

Note. The linear regression of the $\delta^{66}\text{SO}_2/^{64}\text{SO}_2$ values vs. the working standard should be less than **0.06 ‰/V**.

Chapter 8

Technical Information

8.1 Mechanics

Note. This chapter is intended for use by trained Thermo Electron (Bremen) personnel only. Thermo Electron (Bremen) discourages use by and denies liability for the consequences of use by other than Thermo Electron (Bremen) personnel.

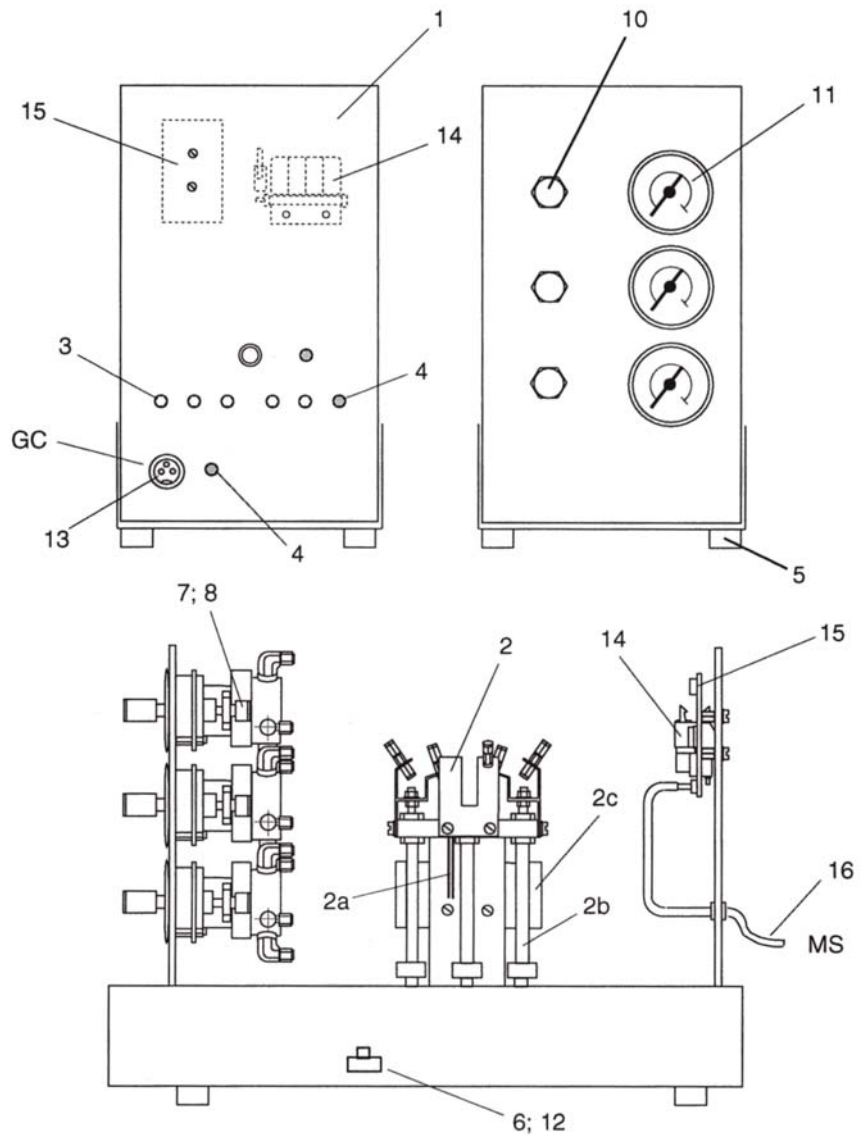


Figure 8-1. ConFlo III Interface (Part No. 113 7200)

Table 8-1. Parts of ConFlo III Interface

Pos. No.	Quantity	Designation	Part No.
1	1	housing	113 7190
2	1	combi inlet	113 2260
2a	2	glass tube with funnel	113 2270
2b	3	lifting cylinder, 8-25	067 4580
2c	1	valve, 5-2 port	113 2350
3	5	coupling	056 7350
4	2	coupling	067 4651
5	4	rubber plate, GM	050 0520
6	1	T-piece, M 5	052 2330
7	3	coupling, 1/8"	074 3360
8	3	ferrule, 1/8"	067 4790
9	1	installation kit ConFlo III	113 7250
10	3	pressure regulator	067 4880
11	3	manometer, 0 -2.5 bar	052 4391
12	2	gasket, 8 * 5	050 5260
13	1	flange type receptacle, BT 3263	032 3760
14	1	manifold card, 4 station, 10 PO	108 3241
15	1	EA start pcb	204 2350
16	1	connection cable	104 0351
17	2 m	silicon hose, 1.0 * 1.75 NF	101 5830

Pos. 9 and 17 are not shown in Figure 8-1.

8.2 Capillaries and Ferrules

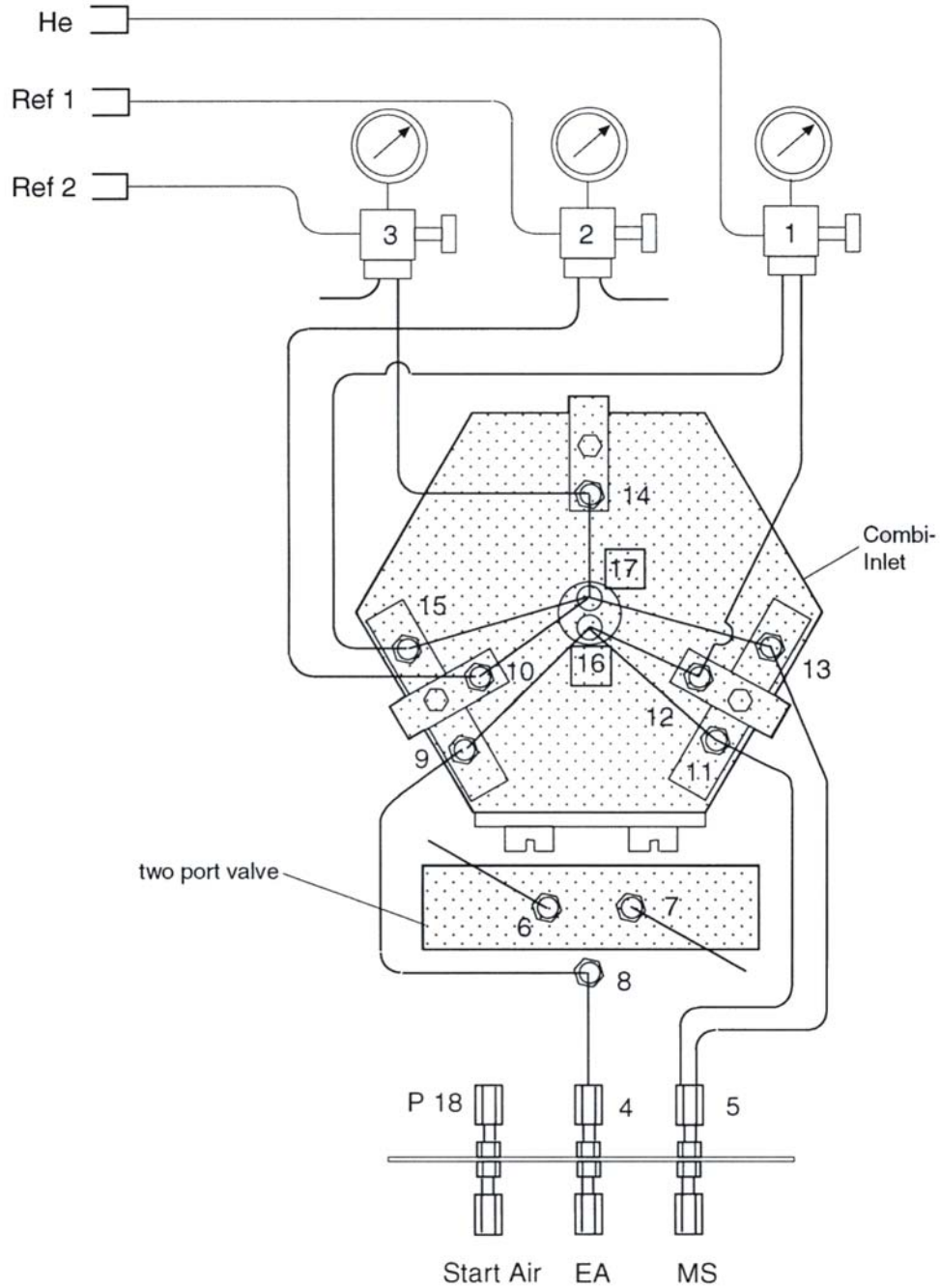


Figure 8-2. Combi Inlet of ConFlo III (Part No. 113 2260)

Table 8-2. Capillaries and Ferrules

Connection	Inner Diameter of <i>Capillary</i>	Length	Part No.	capillary-related <i>Ferrule</i>	Part No.
1 - 12* - 16	0.18 mm	2000 mm	054 2910	GVF2 / 003	106 0170
1 - 15* - 17	0.1 mm	250 mm	104 8990	GVF2 / 003	106 0170
2 - 10* - 17	0.05 mm	310 mm	054 3380	GVF2 / 005	100 4780
3 - 14* - 17	0.05 mm	340 mm	054 3380	GVF2 / 005	100 4780
8 - 9* - 16	0.32 mm	1000 mm	100 4640	GVF / 005	056 6390
4 - 8	1/16"-SS	fitting	060 5470	Swagelock	
5 - 11* - 16	0.1 mm	2000 mm	104 8990	GVF / 003	100 4850
5 - 13* - 17	0.05 mm	1500 mm	054 3380	GVF / 003	100 4850
6	0.32 mm	115 mm	100 4640	GVF / 005	056 6390
7	1/16"-SS	650 mm	060 5470	Swagelock	
2	0.1 mm	200 mm	104 8990	GVF2 / 003	106 0170
3	0.1 mm	200 mm	104 8990	GVF2 / 003	106 0170

* denotes the respective point of support for the capillaries

8.3 Function Schematic

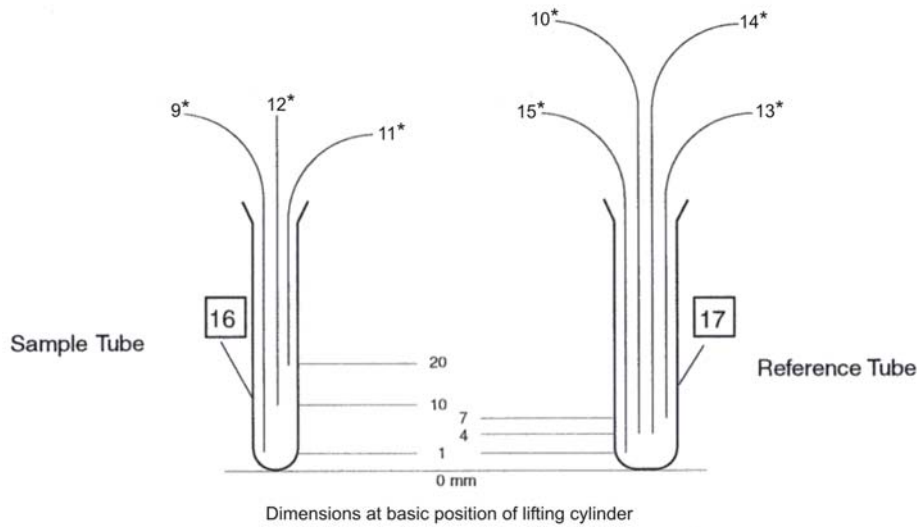


Figure 8-3. Sample Tube and Reference Tube

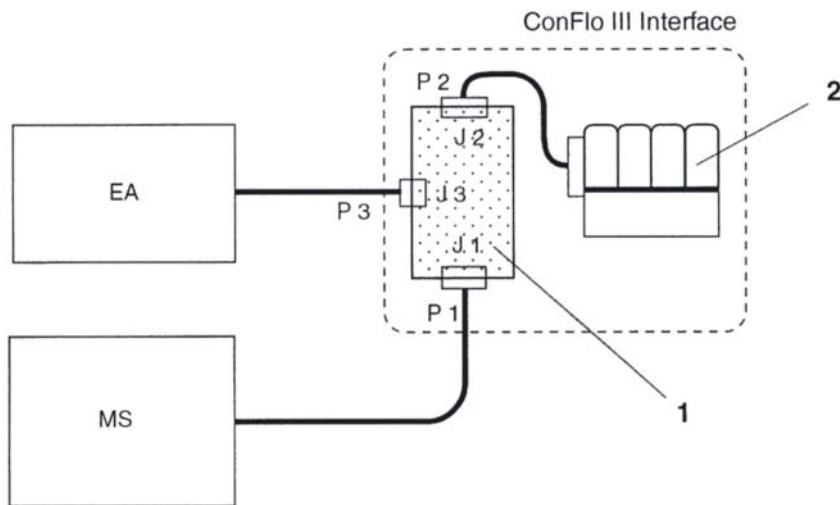


Figure 8-4. Circuit Diagram

Table 8-3. EA Start pcb and Manifold Card in the Circuit Diagram

Pos. No.	Quantity	Designation	Part No.
1	1	EA start pcb	204 2350
2	1	manifold card, 4 station	108 3241

8.4 Connections for Compressed Air

- P 1 - P 18
- P 2 - P 9
- P 3 - P 10
- P 4 - P 5
- P 6 - P 11
- P 7 - P 8

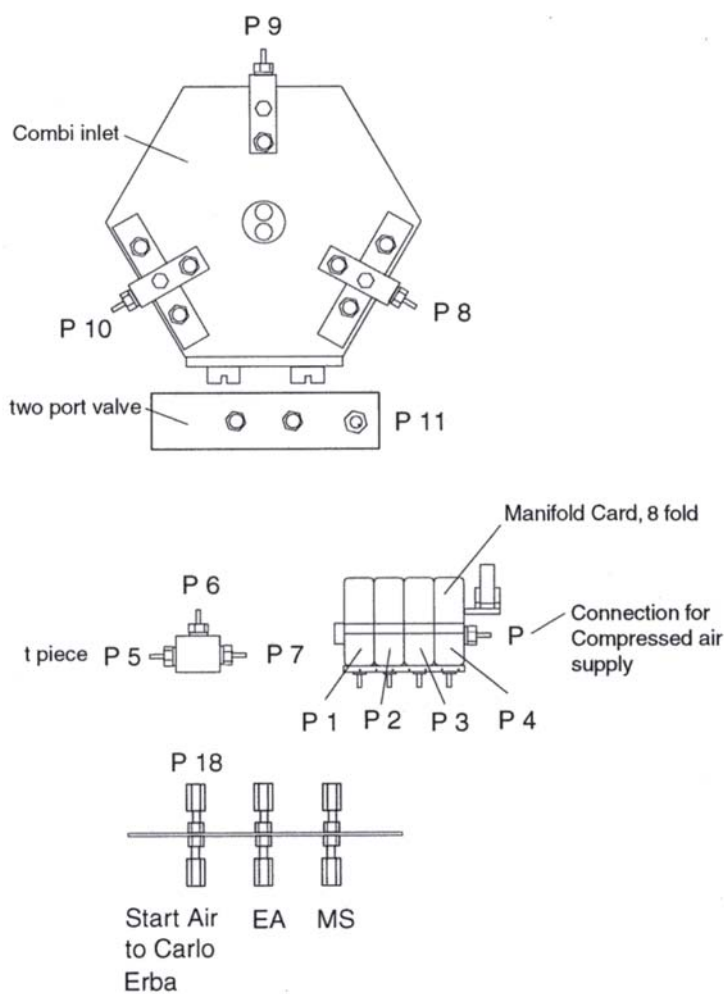


Figure 8-5. Connections for Compressed Air

Index

#

032 3760, 8-3
 049 319, 2-12
 050 0520, 8-3
 050 5260, 8-3
 052 1500, 2-12
 052 2330, 8-3
 052 3210, 2-12
 052 3440, 2-12
 052 3610, 2-12
 052 43 91, 2-12
 052 4391, 8-3
 054 2910, 2-10, 8-5
 054 3380, 2-10, 8-5
 056 6390, 2-11, 8-5
 056 7350, 8-3
 060 5470, 7-4, 8-5
 067 4580, 8-3
 067 4651, 8-3
 067 4790, 2-12, 8-3
 067 4880, 8-3
 074 3360, 2-12, 8-3
 100 4640, 2-10, 8-5
 100 4780, 2-12, 8-5
 100 4850, 2-11, 8-5
 101 5830, 8-3
 104 0351, 8-3
 104 5480, 2-10
 104 8990, 2-10, 8-5
 106 0170, 2-12, 8-5
 108 3241, 8-3, 8-6
 111 8121, 7-4, 7-8
 112 1370, 7-4
 112 1390, 7-4
 113 2260, 8-3, 8-4
 113 2270, 8-3
 113 2350, 8-3
 113 7190, 8-3
 113 7200, 8-2
 113 7250, 8-3
 114 1170, 7-4
 114 1180, 7-4
 114 1210, 7-4
 114 1220, 7-4
 114 1230, 7-4
 115 7100, 7-4
 115 7110, 7-4
 12C16O16O, 5-6
 14N14N, 4-6, 4-7
 15N, 7-2
 15N14N, 4-7
 204 2350, 8-3, 8-6
 204 8580, 7-4
 32S16O16O, 7-12
 32SO₃, 7-3
 34S, 7-3
 34S measurement, 7-12
 34S16O16O, 7-12

34SO₃, 7-3

A

A200S, 3-4, 3-5
 accessories toolbar, 3-8, 3-9, 6-4, 6-5
 accessories window, 3-9, 6-5
 acquisition, 2-5, 4-9, 5-9, 6-3, 7-12, 7-22
 end time, 5-3
 nitrogen, 6-3
 start, 5-3, 7-16
 acquisition end time, 4-3, 5-3, 7-16
 acquisition module, 7-12
 acquisition start, 5-3, 7-16
 acquisition start time, 4-3
 actuator, 2-7
 add configuration, 3-3
 administrative panels button, 3-8, 6-5
 advanced mode, 3-3, 3-4, 3-6
 advanced parameter, 4-5, 5-5
 aerosol spray, 1-2
 air
 atmospheric, 3-11
 compressed, 3-10, 7-9, 8-7
 Al olive, 7-4
 alarm, 7-5
 alcohol, 1-2
 amplification, 7-2
 analysis, 7-8
 carbon, 7-9
 nitrogen, 7-9
 sulfur, 7-2
 analytical condition, 4-2, 5-2, 7-15
 application, 3-12
 Ar, 3-12
 area
 sniffed, 2-4
 AS 128, 3-11
 AS 200, 3-11
 atmospheric air, 3-11
 atmospheric nitrogen, 3-10
 attachment for exhaust tube, 7-4
 autosampler, 3-3, 3-4, 3-5, 3-6, 3-7, 3-10, 3-11, 7-3, 7-9, 7-10
 GC PAL, 3-4, 3-5
 liquid, 3-3, 3-4
 autosampler start signal, 3-8, 6-5

B

back ferrule, 2-12
 background, 3-11, 3-13, 7-10, 7-11
 background value, 3-11, 3-13, 7-10, 7-11
 Bailey, S.A., 7-3
 barium sulfate, 7-3
 basic unit, 8-2
 biological sulfur, 7-2
 board
 driver, 3-10

Finnigan ConFlo III

- bulk sample, 1-4
- button
 - administrative panels, 3-8, 6-5
 - new, 6-7
 - recalibrate, 6-7
 - start, 4-6, 5-6, 7-19
- C**
- cable, 3-10
 - connection, 8-3
 - RS 232, 3-4
- calibration, 7-14
 - jump, 6-4, 6-6, 6-7, 6-8, 6-9
 - mass, 6-3, 7-12
- calibration file, 7-14
- cap nut, 2-12
- capillary, 2-2, 2-4, 2-5, 2-10, 2-11, 3-8, 6-5, 8-4, 8-5
 - carrier gas, 2-4
 - delivering, 2-4
 - dilution, 2-4
 - fixed sample, 2-4
 - fused silica, 3-10, 3-12
 - helium dilution, 3-11, 7-10
 - IRMS, 2-4, 2-5, 2-8
 - Ref 1, 3-8, 6-5
 - Ref 2, 3-8, 6-5
 - reference, 2-4, 2-5
 - sniffer, 2-2, 2-4
 - stainless steel, 3-12
- capillary port, 3-3
- capillary tube, 7-4
- capsule
 - tin, 7-3
- carbon, 1-4, 3-2, 6-2, 6-3, 7-2
 - linearity test, 5-9
 - zero enrichment, 5-2, 5-3, 5-4, 5-5, 5-6, 5-7
- carbon analysis, 7-9
- carbon dioxide, 2-5
- carbon measurement, 5-1, 7-11
- carrier, 7-10
- carrier gas, 2-2, 2-8
- carrier gas capillary, 2-4
- cell, 2-2
 - open split, 2-4, 2-5
- change
 - temperature, 6-4
- change over valve, 7-10
- check, 4-2, 5-2, 7-15
 - leak, 3-10, 3-11, 3-12, 7-9
- chromatogram, 4-3, 4-9, 5-3, 5-6, 5-9, 6-3, 7-16, 7-19, 7-22
- circuit diagram, 8-6
- circuits
 - internal, 1-2
- CO, 2-9, 7-5
- CO₂, 3-8, 3-10, 3-11, 6-5, 6-7, 6-8, 6-9, 7-12, 7-13
- CO₂ configuration, 6-3, 6-4
- CO₂_zero.met, 5-2, 5-9
- collector
 - universal triple, 7-14
- column, 1-4, 3-12, 7-10
 - conditioning, 3-12
 - GC, 3-12, 6-3, 7-9
 - reference 1 - on, 5-2
 - reference 2 - on, 4-2, 7-15
- column conditioning, 3-12
- column temperature, 7-10, 7-11
- COM 1, 3-6
- COM port, 3-4, 3-5, 3-6, 3-7
 - number, 3-7
- COM port number, 3-7
- COM port setting, 3-3, 3-4, 3-5, 3-7
- combi inlet, 2-7, 3-10, 8-3, 8-4
- combustion, 7-2, 7-3
- combustion reactor, 7-2, 7-4
- combustion tube, 7-3
- command
 - properties, 3-6
- communication hardware part, 3-7
- communication parameter, 3-5
- comport, 3-7
- compressed air, 3-10, 7-9, 8-7
- compressed air inlet, 2-8
- compressed air outlet, 2-8
- computer, 6-4
- configuration, 3-3, 3-8, 6-4
 - add, 3-3
 - CO₂, 6-3, 6-4
 - ConFlo, 6-4
 - ConFlo III, 3-10, 7-9
 - cup, 7-14
 - gas, 6-3, 6-4, 6-8, 7-12, 7-13, 7-14
- configurator, 3-3, 3-4
- configurator window, 3-3
- ConFlo configuration, 6-4
- ConFlo II, 2-2
- ConFlo II III interface window, 3-8, 6-5
- ConFlo II set, 3-3
- ConFlo III/III, 3-3
- ConFlo II/III interface, 3-3, 3-9, 6-5
- ConFlo II/III interface window, 3-8, 3-9, 6-5, 6-6
- ConFlo III, 1-2, 1-3, 1-4, 2-6, 2-7, 2-8, 3-2, 3-3, 3-8, 3-10, 3-11, 3-12, 4-2, 4-6, 4-9, 5-2, 5-6, 5-9, 7-4, 7-15, 7-19, 7-22, 8-2, 8-3, 8-4
 - configuration, 3-3
 - installation kit, 8-3
- ConFlo III configuration, 3-3, 3-10, 7-9
- ConFlo III device window, 3-10
- ConFlo III installation, 3-2
- ConFlo III Interface, 2-1, 2-2
- ConFlo III interface, 3-1, 3-8, 4-2, 5-2, 7-5, 7-15, 8-2, 8-3
- ConFlo III set, 3-4
- ConFlo interface, 3-3, 4-2, 5-2, 7-15
- connection, 3-10, 3-12, 7-9, 8-5
 - 1 - 12* - 16, 8-5
 - 1 - 15* - 17, 8-5
 - 2, 8-5
 - 2 - 10* - 17, 8-5
 - 3, 8-5
 - 3 - 14* - 17, 8-5
 - 4 - 8, 8-5
 - 5 - 11* - 16, 8-5
 - 5 - 13* - 17, 8-5

Finnigan ConFlo III

- 6, 8-5
- 7, 8-5
- 8 - 9* - 16, 8-5
- EA/IRMS, 2-2
- connection cable, 8-3
- connection line, 3-12
- connections
 - compressed air, 8-7
- container, 7-3
- continuous flow, 1-4, 2-4
- continuous flow interface, 1-4
- control
 - hardware, 3-7
 - liquid autosampler, 3-5
- convert, 1-4
- copper, 7-3
- copper oxide, 7-8
- coupling, 2-2, 2-12, 8-3
 - 1/8", 8-3
- CuO, 7-3
- cup, 3-11, 3-13, 6-3, 7-11, 7-12, 7-14
 - mass 66, 7-2, 7-10, 7-11
- cup configuration, 7-14
- cycle run time, 7-10

D

- d value, 7-3, 7-12
- d13C/12C, 5-7, 5-9
- d15N, 4-7
- d15N/14N, 4-9
- damage, 1-2
- data, 4-6, 4-9, 5-6, 5-9, 7-19, 7-22
- default setting, 3-7
- default value, 6-8
- delay
 - post jump, 6-8
 - pre jump, 6-8
 - sampling, 7-10
- delivering capillary, 2-4
- Delta IRMS, 7-2
- delta mass spectrometer, 7-14
- detection, 4-9, 5-9, 7-22
- detector
 - H₂, 7-5
- determination
 - on-line, 2-2
 - sulfur, 7-3
- device, 3-5
- devices tab, 3-5
- diagram
 - circuit, 8-6
 - gas flow, 2-9
- diameter
 - inner, 2-4
- dilution, 2-4, 3-8, 6-5
 - helium, 2-10, 2-12, 3-12, 7-10
- dilution capillary, 2-4
- double hole ferrule, 2-12
- driver board, 3-10
- dry out, 2-4
- dual measurement, 6-1, 6-2, 6-3, 6-4

- dynamic flash combustion method, 7-3

E

- EA, 2-10
- EA 1108, 7-9
- EA 1110, 7-3, 7-9
- EA 2500, 7-9
- EA start pcb, 8-3, 8-6
- EA/IRMS connection, 2-2
- Eager software, 3-10, 7-10
- edit menu, 3-4
- editor
 - gas configuration, 7-12
- effluent stream, 2-4
- elbow, 2-12
- elemental analyzer, 1-3, 1-4, 2-2, 2-4, 2-8, 3-1, 3-3, 3-10, 3-11, 3-12, 7-2, 7-3, 7-6, 7-9
 - flash, 7-4
 - manual, 3-10
 - part, 3-12
 - set up, 7-9
- elemental analyzer inlet, 2-8
- elemental analyzer manual, 3-10, 3-12
- elemental analyzer part, 3-12
- elemental analyzer start signal, 3-8, 6-5
- entry, 4-2, 5-2, 7-15
- evaluation tab, 4-4, 5-4, 7-17
- exhaust tube, 7-5, 7-6, 7-9
 - attachment, 7-4
- external surface, 1-2

F

- ferrule, 2-11, 2-12, 8-4, 8-5
 - 1/8", 8-3
 - back, 2-12
 - double hole, 2-12
 - front, 2-12
 - GVF / 003, 8-5
 - GVF / 005, 8-5
 - GVF2 / 003, 8-5
 - GVF2 / 005, 8-5
- field
 - magnetic, 6-3
- file
 - calibration, 7-14
- file browser, 4-2, 5-2, 7-15
- final temperature, 3-12
- fixed sample capillary, 2-4
- flange type receptacle, 8-3
- flash combustion, 7-3
- flash elemental analyzer, 7-4, 7-10
- flow, 2-4, 3-12
 - helium, 3-11, 3-12
 - sample, 2-4
- focus setting, 3-11, 3-13
- front ferrule, 2-12
- front panel, 2-6
- function schematic, 8-6
- funnel, 8-3

furnace
 oxidation, 3-12
 reduction, 1-4, 3-12
fused silica, 2-8, 2-10, 2-11
fused silica capillary, 3-10, 3-12

G

gas, 1-4, 2-4, 2-5, 2-9, 6-3, 7-5, 7-12
 reference, 2-12, 4-2, 5-2, 6-7, 7-15
 standard, 3-12
 toxic, 2-9
gas configuration, 6-3, 6-4, 6-8, 7-12, 7-13, 7-14
gas configuration editor, 7-12
gas connection, 2-8
gas flow, 2-2
gas flow diagram, 2-9
gas introduction, 2-2
gas isotope mass spectrometer, 1-4
gas pressure, 3-10, 3-11, 7-9
gas supply, 3-10
gas tank, 3-10
gaschromatographic, 1-4
gasket, 8-3
GC column, 3-12, 6-3, 7-9
 SO₂, 7-6
 stainless steel, 7-9, 7-10
 Teflon, 7-4
 teflon, 7-9, 7-11
GC PAL autosampler, 3-4, 3-5
GC/ConFlo III installation menu, 3-2
GEM installation menu, 3-2
glass tube, 8-3
guide
 user, 1-3

H

H₂, 7-5
H₂ detector, 7-5
H₂O, 7-10, 7-11
hardware, 3-3
hardware control, 3-7
hazardous, 7-5
health, 7-5
heat, 1-2
heater
 ion source, 3-12, 7-10
 needle valve, 3-12
heating foil, 7-4
helium, 2-2, 2-4, 3-10, 3-11, 7-3
helium dilution, 2-4, 2-10, 2-12, 3-12, 7-10
helium dilution capillary, 3-11, 7-10
helium flow, 3-11, 3-12
helium inlet, 2-8
helium pressure gauge, 2-6
helium pressure regulator, 2-6
helium purge, 2-10, 2-12
helium stream, 7-3
high temperature conversion, 1-4
high vacuum, 3-11

high voltage setting, 6-7
hole
 single, 2-11
hose
 silicon, 8-3
housing, 8-3
hydrogen, 1-4, 3-2

I

ignition, 1-2
information
 technical, 2-7, 8-1
inlet
 combi, 2-7, 3-10, 8-4
 compressed air, 2-8
 elemental analyzer, 2-8
 helium, 2-8
 reference, 2-2
 reference gas, 3-8, 6-5
 reference gas 1, 2-8
 reference gas 2, 2-8
 sample gas, 2-2
inlet system, 1-4
inner diameter, 2-4
installation, 3-4, 6-2
 ConFlo III, 3-2
instrument control, 3-8, 6-4
instrument control window, 6-6
instrument tab, 4-2, 5-2, 7-15
interface
 ConFlo, 4-2, 5-2, 7-15
 ConFlo III, 4-2, 5-2, 7-5, 7-15
 RS 232, 3-3
interface type, 3-7
internal circuits, 1-2
introduction
 gas, 2-2
ion intensity, 3-13, 4-6, 5-6, 7-19
ion source, 1-4, 2-4, 3-11, 7-10
ion source heater, 3-12, 7-10
IRMS, 1-3, 1-4, 2-2, 2-10, 3-10, 3-12, 3-13, 4-2, 5-2, 6-2,
7-9, 7-12, 7-14, 7-15
 Delta, 7-2
IRMS capillary, 2-4, 2-5, 2-8
IRMS connection cable, 2-8
Isodat 2.0, 1-3, 3-3, 3-4, 3-10
 setup, 3-10
Isodat 2.0 setup, 3-10
Isodat installation menu, 3-2
Isodat mode, 3-2
Isodat old, 3-2
isotope, 7-2
 ratio, 1-4
isotope ratio, 1-3, 1-4, 3-2, 7-2, 7-3
isotope ratio analysis, 2-2
 on-line, 1-4
isotope ratio mass spectrometer, 1-3, 1-4, 2-2
isotope ratio measurement, 7-2

Finnigan ConFlo III

J

jump
 magnet, 6-3
 jump calibration, 6-4, 6-6, 6-7, 6-8, 6-9

K

kit
 installation ConFlo III, 8-3
 sulfur measurement, 7-4

L

leak, 3-13
 leak check, 3-10, 3-11, 3-12, 7-9
 left side view, 2-7
 liability, 8-2
 lifting cylinder, 8-3
 line
 connection, 3-12
 linear regression, 4-9, 5-9, 7-23
 linearity test, 4-2, 5-2, 7-15
 carbon, 5-9
 nitrogen, 4-9
 sulfur, 7-22, 7-23
 liquid autosampler, 3-3, 3-4
 liquid autosampler control, 3-5
 list
 time events, 4-2, 4-4, 5-2, 5-4, 5-6, 7-15, 7-17

M

magnet field value, 7-12
 magnet jump, 6-3
 magnet setting, 6-7
 magnetic field, 6-3
 main power supply, 3-10
 main valve, 3-10
 manifold card, 2-7, 8-3, 8-6
 manometer, 2-11, 2-12, 3-12, 7-7, 8-3
 part, 2-12
 rear side, 7-7
 manometer part, 2-12
 manual, 1-3
 changes, vi
 elemental analyzer, 3-12
 manual operation, 3-8
 MAS 200, 3-11
 mass, 3-11, 3-13, 7-11, 7-12, 7-14
 nitrogen, 6-3
 predefined, 7-12
 mass 28, 4-6
 mass 40, 3-12
 mass 44, 5-6, 6-3, 7-14
 mass 45, 6-3, 7-14
 mass 46, 6-3, 7-14
 mass 64, 7-12, 7-14, 7-19
 mass 66, 7-2, 7-12, 7-14
 mass calibration, 6-3, 7-12
 mass spectrometer, 2-2, 2-5

delta, 7-14
 gas isotope, 1-4
 material
 organic, 1-4, 3-12
 matter
 organic, 3-2
 Max/Min setting, 6-7
 measurement, 1-1, 3-12, 4-6, 4-9, 5-6, 5-9, 7-14
 34S, 7-12
 carbon, 5-1, 7-11
 dual, 6-2, 6-3, 6-4
 isotope ratio, 7-2
 nitrogen, 4-1, 7-11
 single element, 6-4
 sulfur, 7-1, 7-2, 7-3, 7-4, 7-6, 7-9, 7-12
 mechanics, 8-2
 menu
 edit, 3-4
 GC/ConFlo III installation, 3-2
 GEM installation, 3-2
 Isodat installation, 3-2
 method, 4-2, 4-9, 5-2, 5-9, 6-3, 7-15, 7-22
 dynamic flash combustion, 7-3
 mixture, 7-2
 organic and inorganic compound, 7-2
 mode
 advanced, 3-3, 3-4, 3-6
 Isodat, 3-2
 normal, 3-6
 module
 acquisition, 7-12

N

N₂, 3-10, 3-11, 6-3, 6-4, 6-7, 6-8, 6-9
 N₂_zero.met, 4-2
 needle valve, 3-10, 3-11, 7-10
 needle valve heater, 3-12
 new button, 6-7
 nitrogen, 1-4, 2-5, 3-2, 4-2, 4-3, 4-4, 4-5, 4-6, 6-2, 6-3, 7-2, 7-11
 atmospheric, 3-10
 linearity test, 4-9
 zero enrichment, 4-2, 4-7
 nitrogen acquisition, 6-3
 nitrogen analysis, 7-9
 nitrogen mass, 6-3
 nitrogen measurement, 4-1, 7-11
 nitrogen peak, 6-3
 nitrogen trace, 7-3
 normal mode, 3-6

O

O₂, 3-10, 3-11, 7-3, 7-9
 O₂ pressure, 7-3
 off-time, 4-3, 5-3, 7-16
 off-value, 4-4, 5-4, 7-17
 old Isodat, 3-2
 olive
 Al, 7-4

on-line determination, 2-2
 on-line isotope ratio analysis, 1-4
 on-off, 4-2, 5-2, 7-15
 on-off entry, 4-2, 5-2, 7-15
 on-off pulse, 4-3, 5-3, 5-6, 7-16
 open split arrangement, 2-2, 2-4
 open split cell, 2-2, 2-4, 2-5, 3-11
 open split principle, 2-2
 operating signal, 3-4
 operation
 manual, 3-8
 valve, 2-4
 operational safety, 7-5
 organic material, 1-4, 3-12
 organic matter, 3-2
 organic solvent, 1-2
 outlet
 compressed air, 2-8
 oxidation, 1-4
 oxidation furnace, 3-12
 oxidation reactor, 3-10
 oxygen, 1-4, 3-2, 7-3, 7-10

P

P 1, 8-7
 P 10, 8-7
 P 11, 8-7
 P 18, 8-7
 P 2, 8-7
 P 3, 8-7
 P 4, 8-7
 P 5, 8-7
 P 6, 8-7
 P 7, 8-7
 P 8, 8-7
 P 9, 8-7
 packing, 7-8
 reactor, 7-8
 paint thinner, 1-2
 panel
 front, 2-6
 rear, 2-6, 2-8, 2-9, 3-10
 parameter
 advanced, 4-5, 5-5
 communication, 3-5
 peak, 4-7, 4-9, 5-6, 5-7, 5-9, 6-9, 7-20, 7-22
 nitrogen, 6-3
 reference gas, 4-4, 5-4, 6-3, 7-17
 sample, 6-3
 peak center, 6-4, 6-9
 peak detection tab, 4-5, 5-5, 7-18
 piston, 2-4, 2-5, 2-7, 3-10
 plate
 rubber, 8-3
 pneumatic valve, 2-4
 point of support, 8-5
 port
 capillary, 3-3
 COM, 3-4, 3-6, 3-7
 reference, 4-2, 5-2, 6-8, 7-15
 reference gas, 4-2, 5-2, 7-15

 source, 3-3
 post jump delay, 6-8
 power supply, 7-4
 pre jump delay, 6-8
 precaution, 1-2
 precautions, 1-2
 precision, 2-2
 predefined mass, 7-12
 prediluter, 2-8
 pressure, 3-10, 3-12
 gas, 3-10, 3-11, 7-9
 O₂, 7-3
 reference gas, 4-9, 5-9, 7-22
 pressure gauge, 2-7
 reference gas 1, 2-6
 reference gas 2, 2-6
 pressure regulator, 4-6, 5-6, 7-19, 8-3
 reference gas 1, 2-6
 reference gas 2, 2-6
 printout tab, 4-6, 5-6, 7-19
 properties command, 3-6
 pulse
 on-off, 4-3, 5-3, 5-6, 7-16
 reference gas, 4-3, 5-3, 7-16
 purge, 3-11, 7-9
 helium, 2-10, 2-12
 reference, 2-12

Q

quality, 3-10

R

radiator, 1-2
 range, 7-10, 7-11
 ratio, 7-12
 ratio group, 7-12, 7-13
 ratio groups window, 7-13
 reactor, 7-3, 7-6, 7-8, 7-9, 7-10, 7-11
 combustion, 7-2, 7-4
 filling, 7-8
 oxidation, 3-10
 ready for use, 7-8
 reduction, 3-10
 SO₂, 7-9
 reactor filling, 7-8
 reactor packing, 7-8
 reactor temperature, 7-11
 ready for use, 7-4
 ready for use reactor, 7-8
 rear panel, 2-6, 2-8, 2-9, 3-10
 recalibrate button, 6-7
 reducing valve, 3-10
 reduction furnace, 1-4, 3-12
 reduction reactor, 3-10
 Ref 1, 3-10
 Ref 1 capillary, 3-8, 6-5
 Ref 2, 3-10
 Ref 2 capillary, 3-8, 6-5
 Ref. 1, 2-10

Finnigan ConFlo III

- Ref. 2, 2-10
 - ref. time, 5-4
 - reference, 2-2, 2-5, 7-10
 - reference 1, 3-8, 5-2, 6-5
 - reference 1 - on column, 5-2
 - reference 2, 3-8, 4-2, 6-5, 7-15
 - reference 2 - on column, 4-2, 7-15
 - reference capillary, 2-4, 2-5
 - reference gas, 2-9, 2-12, 4-2, 5-2, 6-7, 7-15
 - reference gas 1 inlet, 2-8
 - reference gas 1 pressure gauge, 2-6
 - reference gas 1 pressure regulator, 2-6
 - reference gas 2 inlet, 2-8
 - reference gas 2 pressure gauge, 2-6
 - reference gas 2 pressure regulator, 2-6
 - reference gas inlet, 3-8, 6-5
 - reference gas peak, 4-4, 5-4, 6-3, 7-17
 - reference gas port, 4-2, 5-2, 7-15
 - reference gas pressure, 4-9, 5-9, 7-22
 - reference gas pulse, 4-3, 5-3, 7-16
 - reference inlet, 2-2
 - reference port, 4-2, 5-2, 6-8, 7-15
 - reference purge, 2-12
 - reference section, 2-2, 2-3, 2-4, 2-10, 2-11
 - reference time, 4-4, 7-17
 - reference tube, 8-6
 - regulator
 - pressure, 4-6, 5-6, 7-19, 8-3
 - resistor, 3-11, 3-13
 - result grid, 4-9, 5-7, 5-9, 7-20, 7-23
 - right side view, 2-8
 - RS 232 cable, 3-4
 - RS 232 interface, 3-3
 - rubber plate, 8-3
 - run, 6-4
 - single, 5-9, 7-22
- ## S
- safety
 - operational, 7-5
 - sample, 1-4, 2-2, 2-4, 3-12, 4-2, 5-2, 6-2, 6-3, 7-3, 7-15
 - sample flow, 2-4
 - sample gas inlet, 2-2
 - sample peak, 6-3
 - sample preparation, 1-3, 1-4
 - sample section, 2-2, 2-3, 2-4, 2-10, 2-11
 - sample size, 7-2
 - sample stream, 2-2
 - sample tube, 8-6
 - sampling delay, 7-10
 - script, 3-4
 - section
 - reference, 2-3, 2-4, 2-10, 2-11
 - sample, 2-3, 2-4, 2-10, 2-11
 - sensitivity, 3-11, 3-13, 7-10, 7-11
 - separation, 6-3
 - sequence, 4-6, 5-6
 - set
 - ConFlo II, 3-3
 - ConFlo III, 3-4
 - set up
 - elemental analyzer, 7-9
 - setting, 3-3, 3-5, 3-7, 3-10
 - COM port, 3-3, 3-4, 3-5, 3-7
 - default, 3-7
 - focus, 3-11, 3-13
 - high voltage, 6-7
 - magnet, 6-7
 - Max/Min, 6-7
 - side view, 2-7
 - signal
 - operating, 3-4
 - start, 3-10
 - signal height, 6-9
 - silicon hose, 8-3
 - single element measurement, 6-4
 - single hole, 2-11
 - single run, 4-9, 5-9, 7-22
 - size
 - sample, 7-2
 - Smith, J.W., 7-3
 - sniffed area, 2-4
 - sniffer capillary, 2-2, 2-4
 - sniffer point, 2-4
 - sniffing position, 2-5
 - SO₂, 2-9, 7-3, 7-5, 7-7, 7-9, 7-12, 7-13, 7-14
 - condensation, 7-7
 - SO₂ condensation, 7-7
 - SO₂ GC column, 7-6
 - SO₂ reactor, 7-9
 - SO₂_zero.met, 7-15, 7-22
 - SO₃, 7-3
 - software, 1-3, 3-2, 3-4, 3-10
 - Eager, 3-10
 - solvent
 - organic, 1-2
 - source, 3-11, 6-7
 - source port, 3-3
 - spare part, 2-10, 2-11
 - spray, 1-2
 - stainless steel, 3-10
 - stainless steel capillary, 3-12
 - stainless steel GC column, 7-9, 7-10
 - stainless steel tubing, 7-6
 - standard, 2-5
 - standard deviation, 4-7, 5-7, 7-20
 - standard gas, 2-9, 3-12
 - standard on/off test, 4-2, 7-15
 - start button, 4-6, 5-6, 7-19
 - start cable, 2-8
 - start signal, 3-10
 - autosampler, 3-8, 6-5
 - elemental analyzer, 3-8, 6-5
 - stream
 - effluent, 2-4
 - sample, 2-2
 - sulfur, 1-4, 3-2, 7-2
 - biological, 7-2
 - linearity test, 7-22, 7-23
 - zero enrichment, 7-15
 - sulfur analysis, 7-2
 - sulfur determination, 7-3
 - sulfur dioxide, 7-3, 7-5

sulfur measurement, 7-1, 7-2, 7-3, 7-4, 7-6, 7-9, 7-12
sulfur measurement kit, 7-4
sulfur zero enrichment, 7-15, 7-16, 7-17, 7-18, 7-19, 7-20
sunlight, 1-2
supply
 gas, 3-10
 main power, 3-10
 power, 7-4
support
 point of, 8-5
surface
 external, 1-2
Swagelock, 8-5
system, 6-3

T

tab
 devices, 3-5
 evaluation, 4-4, 5-4, 7-17
 instrument, 4-2, 5-2, 7-15
 peak detection, 4-5, 5-5, 7-18
 printout, 4-6, 5-6, 7-19
 time events, 4-3, 5-3, 7-16
TC/EA, 1-4
technical information, 2-7, 8-1
teflon, 2-12, 3-10
Teflon GC column, 7-4
teflon GC column, 7-9, 7-11
Teflon tube, 7-4
teflon tube, 7-11
teflon tubing, 3-10, 7-6, 7-9
temperature, 7-7, 7-10
 column, 7-11
 final, 3-12
 reactor, 7-11
temperature change, 6-4
template, 7-13
test
 linearity, 4-2, 5-2, 7-15
 standard on/off, 4-2, 7-15
Thermo Electron, 1-4, 2-9, 7-4, 7-8
Thermo Electron (Bremen), 7-2, 8-2
time, 6-3, 6-4, 7-8
 acquisition end, 4-3, 7-16
 acquisition start, 4-3
 cycle run, 7-10
 off, 4-3, 5-3, 7-16
 ref., 5-4
 reference, 4-4, 7-17
time events list, 4-2, 4-4, 5-2, 5-4, 5-6, 7-15, 7-17
time events tab, 4-3, 5-3, 7-16
tin capsule, 7-3
toolbar
 accessories, 3-8, 3-9, 6-4, 6-5
toxic gas, 2-9
t-piece, 2-10, 8-3
trap
 water, 3-10
tube
 capillary, 7-4
 combustion, 7-3

 exhaust, 7-5, 7-6, 7-9
 reference, 8-6
 sample, 8-6
 Teflon, 7-4
 teflon, 7-11
tube in tube, 2-2
tubing
 stainless steel, 7-6
 teflon, 3-10, 7-6, 7-9
tungsten oxide, 7-8
two port valve, 2-8

U

union, 7-4
universal triple collector, 7-14
user guide, 1-3

V

vacuum
 high, 3-11
value
 background, 3-11, 3-13, 7-10, 7-11
 d, 7-12
 default, 6-8
 magnet field, 7-12
 off, 4-4, 5-4, 7-17
valve, 3-10, 3-12
 5-2 port, 8-3
 change over, 7-10
 main, 3-10
 needle, 3-10, 3-11, 7-10
 pneumatic, 2-4
 reducing, 3-10
 two port, 2-8
valve operation, 2-4
vent, 2-10
ventilation, 7-5, 7-7, 7-9
view
 left side, 2-7
 right side, 2-8
 side, 2-7

W

water, 3-12
water trace, 7-3
water trap, 3-10, 7-9
window
 accessories, 3-9, 6-5
 configurator, 3-3
 ConFlo II III interface, 3-8, 6-5
 ConFlo II/III interface, 3-8, 3-9, 6-5, 6-6
 ConFlo III device, 3-10
 instrument control, 6-6
 ratio groups, 7-13
working area, 7-5
working standard, 4-9, 5-9, 7-23

Finnigan ConFlo III

Z

Zero Enrichment, 4-3

zero enrichment, 4-2, 4-3, 4-4, 4-5, 4-6, 4-9, 5-2, 5-9,
7-15, 7-22

nitrogen, 4-2

sulfur, 7-15, 7-16, 7-17, 7-18, 7-19, 7-20

